biotechrabbit

Product Catalog 2024–2025







biotechrabbit GmbH Volmerstr.9 12489 Berlin, Germany info@biotechrabbit.com www.biotechrabbit.com Office: +493055578210 Fax: +4930555782199

Managing Director: Dr. Bernd Haase Commercial Register: Charlottenburg HRB 232169 B VAT No/USt-Id: DE 279969868

Bank account Deutsche Bank Berlin BIC/SWIFT: DEUTDEDBBER IBAN: DE22 1007 0024 0062 8891 02



About biotechrabbit

Scientists around the world are working to leap ahead of diseases and to improve our lives with innovation. Aside from brilliant minds and relentless passion, the success of science depends on the quality of the materials used.

biotechrabbit is determined to offer the best products and services to those who lead the way in research. The valued relationships with our partners and customers drive us to exceed current limitations with flexibility, innovation and customized solutions to match their specific requirements. Each member of our team of highly engaged scientists, experienced managers and talented business developers aims to facilitate our partners and customers to leap and lead progress in the life science.

Our way of doing business combines the passion and pure curiosity of excellent researchers with the agile spirit of true entrepreneurs.

biotechrabbit leap and lead

Products made in Germany

- Diagnostic-grade quality
- PCR enzymes and mixes
- Nucleic acid purification
- Monoclonal Antibodies
- Cell-free protein synthesis

Services

- · OEM, custom and bulk
- Lyophilization services

Contact Us

Phone +49 30 555 78 21 10

General contact info@biotechrabbit.com

Ordering order@biotechrabbit.com

Technical Support support@biotechrabbit.com

Sales sales@biotechrabbit.com Fax +49 30 555 78 21 99

OEM and Services oem@biotechrabbit.com

Collaboration and Partnership partnership@biotechrabbit.com

Career Opportunities hr@biotechrabbit.com

Product Overview

Custom Solutions and Diagnostic Services	1
About biotechrabbit	2
Lyophilization	4
Custom Protein Manufacturing	6
Antibody Services	7
Standard PCR Taq DNA Polymerase YourTaq [™] PCR Master Mix YourTaq [™] PCR Master Mix Direct-Load AmPelify [™] PCR Master Mix Green PCR Master Mix Direct-Load	10 11 12 12 13 13
Hot-Start PCR (Antibody-mediated) UPstart [™] Taq Antibody UPstart [™] Plus Taq Antibody YourTaq [™] Hot-Start DNA Polymerase Hot-Start Taq DNA Polymerase YourTaq [™] Hot-Start PCR Master Mix Hot-Start PCR Master Mix YourTaq [™] Hot-Start PCR Master Mix Direct-Load Green Hot-Start PCR Master Mix Direct-Load	15 16 17 17 18 18 18 19 19
Hot-Start PCR (Aptamer-mediated) ApStarTaq [™] Hot-Start DNA Polymerase ApStarTaq [™] Hot-Start PCR Master Mix	<mark>20</mark> 21 21
Multiplex Standard PCR Multiplex Hot-Start PCR Master Mix	<mark>22</mark> 22
Reverse Transcription and RT-PCR AnyScript [™] Reverse Transcriptase AllScript [™] Reverse Transcriptase RevertUP [™] II Reverse Transcriptase MMuLV Reverse Transcriptase 1-Step RT-PCR Master Mix cDNA Synthesis Kit	23 24 25 25 26 26
Quantitative Multiplex Real-Time PCRCAPITAL™ qPCR Probe Master MixCAPITAL™ 1-Step qRT-PCR Probe Master MixCAPITAL™ qPCR Probe Master Mix PlusCAPITAL™ 1-Step qRT-PCR Probe Master Mix PlusCAPITAL™ qPCR Green Master MixCAPITAL™ 1-Step qRT-PCR Green Master MixStar™ qPCR Probe Master MixUniFi™ 1-Step qRT-PCR Probe Master Mix	28 29 30 30 31 31 32 32
LYO-ready Master Mix LYO-ready CAPITAL [™] qPCR Probe Mix LYO-ready CAPITAL [™] 1-Step qRT-PCR Probe Mix LYO-ready CAPITAL [™] qPCR Probe Mix Plus LYO-ready CAPITAL [™] 1-Step qRT-PCR Probe Mix Plus	<mark>34</mark> 35 35 36 36



Product Overview

Lyophilized Polymerases and Master Mixes	38
Hot-Start Taq DNA Polymerase, lyophilized	38
YourTaq [™] Hot-Start PCR Master Mix, lyophilized	39
Hot-Start PCR Master Mix, lyophilized	39
CAPITAL [™] qPCR Probe Master Mix, lyophilized	40
CAPITAL [™] 1-Step qRT-PCR Probe Master Mix, lyophilized	40
Long-Range PCR, High-Fidelity and NGS	42
Long Range DNA Polymerase	43
Long Range PCR Master Mix	43
Pfu DNA Polymerase	44
Pfu PCR Master Mix	44
T4 DNA Ligase Rapid	45
ExactiFi [™] Hot-Start DNA Polymerase	45
A-Tailing Kit	46
DNA Blunting Kit	46
Enzymes and Proteins for Molecular Biology	48
LYO-ready T6 UvsX DNA Recombinase	49
LYO-ready T4 UvsY Protein	49
LYO-ready Exonuclease III	50
LYO-ready RB69 Gene 32 Protein	50
LYO-ready T4 Gene 32 Protein	51
RNase Inhibitor	51
Heat Labile Uracil-DNA Glycosylase	52
Proteinase K	52
T7 RNA Polymerase	53
phi29 DNA Polymerase	53
Nucleotides, Additives and Accessories	54
Nucleotides, Additives and Accessories dNTP Sets and Mixes	<mark>54</mark> 54
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂	<mark>54</mark> 54 55
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer	<mark>54</mark> 54 55 55
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water	<mark>54</mark> 54 55 55 55
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification	54 54 55 55 55
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup	54 54 55 55 55 55
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit	54 54 55 55 55 56 58
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit	54 54 55 55 55 56 58 58
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation	54 55 55 55 56 58 58
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl2 PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit	54 55 55 55 56 58 58 58 58
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation	54 55 55 55 56 58 58 58
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation GenUP [™] Bacterial gDNA Kit	54 55 55 55 56 58 58 58 59 60
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation GenUP [™] Bacterial gDNA Kit GenUP [™] gDNA Kit	54 55 55 55 56 58 58 58 59 60 60
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit Total DNA Isolation GenUP™ Bacterial gDNA Kit GenUP™ Blood DNA Kit	54 55 55 55 56 58 58 58 59 60 60 60 61
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit Total DNA Isolation GenUP™ Bacterial gDNA Kit GenUP™ Blood DNA Kit GenUP™ Plant DNA Kit	54 55 55 55 56 58 58 58 59 60 60 61 61
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation GenUP [™] Blacterial gDNA Kit GenUP [™] Blood DNA Kit GenUP [™] Blood DNA Kit Total RNA Isolation	54 55 55 56 58 58 58 59 60 60 61 61
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation GenUP [™] Blacterial gDNA Kit GenUP [™] Blood DNA Kit GenUP [™] Plant DNA Kit Total RNA Isolation GenUP [™] Total RNA Kit	 54 55 55 56 58 58 59 60 60 61 61 62
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit Total DNA Isolation GenUP™ Bacterial gDNA Kit GenUP™ Blood DNA Kit GenUP™ Blood DNA Kit GenUP™ Plant DNA Kit Total RNA Isolation GenUP™ Total RNA Kit GenUP™ Micro RNA Kit	 54 54 55 55 56 58 58 59 60 60 61 61 62 62 62
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation GenUP [™] Bacterial gDNA Kit GenUP [™] Blood DNA Kit GenUP [™] Blood DNA Kit GenUP [™] Plant DNA Kit GenUP [™] Plant DNA Kit GenUP [™] Plant DNA Kit GenUP [™] Micro RNA Kit GenUP [™] Micro RNA Kit GenUP [™] Blood RNA Kit	 54 55 55 56 58 58 59 60 60 61 61 62 62 63
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit Total DNA Isolation GenUP™ Bacterial gDNA Kit GenUP™ Blood DNA Kit GenUP™ Blood DNA Kit GenUP™ Plant DNA Kit GenUP™ Plant DNA Kit GenUP™ Plant AKit GenUP™ Micro RNA Kit GenUP™ Blood RNA Kit GenUP™ Plant RNA Kit	 54 54 55 55 56 58 58 59 60 60 61 61 62 63 63
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP [™] PCR/Gel Cleanup Kit GenUP [™] Exo SAP Kit Plasmid Isolation GenUP [™] Plasmid Kit Total DNA Isolation GenUP [™] Bacterial gDNA Kit GenUP [™] Blood DNA Kit GenUP [™] Blood DNA Kit GenUP [™] Total RNA Kit GenUP [™] Total RNA Kit GenUP [™] Micro RNA Kit GenUP [™] Blood RNA Kit GenUP [™] Plant RNA Kit Kirus DNA and RNA Isolation	 54 55 55 56 58 58 59 60 60 61 61 62 63 63
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit Total DNA Isolation GenUP™ Bacterial gDNA Kit GenUP™ Blood DNA Kit GenUP™ Plant DNA Kit Total RNA Isolation GenUP™ Diant RNA Kit GenUP™ Micro RNA Kit GenUP™ Blood RNA Kit GenUP™ Plant RNA Kit Virus DNA and RNA Isolation GenUP™ BS Virus RNA Kit	 54 55 55 56 58 58 59 60 60 61 61 62 63 63 64
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP™ PCR/Gel Cleanup Kit GenUP™ Exo SAP Kit Plasmid Isolation GenUP™ Plasmid Kit Total DNA Isolation GenUP™ Bacterial gDNA Kit GenUP™ Blood DNA Kit GenUP™ Plant DNA Kit Total RNA Isolation GenUP™ Plant DNA Kit GenUP™ Plant RNA Kit GenUP™ Micro RNA Kit GenUP™ Blood RNA Kit GenUP™ Plant RNA Kit GenUP™ Plant RNA Kit GenUP™ Blood RNA Kit	 54 55 55 56 58 58 59 60 61 61 62 63 63 64 64
Nucleotides, Additives and Accessories dNTP Sets and Mixes MgCl ₂ PCR Enhancer PCR Grade Water Nucleic Acid Purification DNA Cleanup GenUP ^{IM} PCR/Gel Cleanup Kit GenUP ^{IM} Exo SAP Kit Plasmid Isolation GenUP ^{IM} Plasmid Kit Total DNA Isolation GenUP ^{IM} Blood DNA Kit GenUP ^{IM} Blood DNA Kit GenUP ^{IM} Plant DNA Kit GenUP ^{IM} Plant DNA Kit GenUP ^{IM} Plant DNA Kit GenUP ^{IM} Micro RNA Kit GenUP ^{IM} Blood RNA Kit GenUP ^{IM} Blood RNA Kit GenUP ^{IM} Blood RNA Kit GenUP ^{IM} Plant RNA Kit GenUP ^{IM} Blood RNA Kit	 54 54 55 55 56 58 58 59 60 61 61 62 63 63 64 64

Product Overview

Electrophoresis	66
DNA Electrophoresis Ladders and Loading Dye	66
Protein Electrophoresis Ladders	67
Cell-free Protein Synthesis	69
E. coli Protein Synthesis Kits	
RTS™ 100 <i>E. coli</i> HY Kit	70
RTS™ 500 ProteoMaster <i>E. coli</i> HY Kit	70
RTS™ 9000 <i>E. coli</i> HY Kit	71
RTS™ 100/500 <i>E. coli</i> Disulfide Kits	71
Wheat Germ Protein Synthesis Kits	
RTS [™] 100 Wheat Germ Kit	72
RTS [™] 100/500 Wheat Germ CECF Kits	72
Insect Protein Synthesis Kit	
RTS [™] 100 Insect Membrane Kit	73
RTS Supplements	
RTS [™] Amino Acid Sampler	73
RTS™ DnaK Supplement	74
RTS [™] GroE Supplement	74
Linear Templates and Vectors for Cell-Free Protein Synthesis	75
Linear Templates for Cell-free Protein Synthesis	
RTS [™] Linear Template Kit Plus	76
RTS [™] Wheat Germ LinTempGen Set, His ₆ -tag	76
RTS™ Linear Template Fab Kit	77
RTS™ 100 <i>E. coli</i> LinTempGen Set, His-tag	77
Vectors for Cell-free Protein Synthesis	
RTS™pIX4.0 Insect Vector	78
RTS™pIX3.0 Vector	78
RTS [™] pIVEX Wheat Germ His6-tag Vector Set	79
RTS [™] pIVEX <i>E. coli</i> His-tag, 2nd Gen. Vector Set	79
High Affinity His-tag Binding	80
Tris-NTA Amine	80
Tris-NTA Biotin	80



Custom Solutions and Diagnostic Services

biotechrabbit innovates, develops and manufactures products with superior performance for diagnostics, life science research and applied markets. We specialize in the production of enzymes and kits for molecular biology and provide high-quality products with top-class availability.

biotechrabbit products are available in bulk or custom formulations, matching each customer's individual needs.

- Enzymes for molecular applications
- Contract manufacturing
- · Lyophilization services
- OEM and Private Label

biotechrabbit's collaborative approach ensures reliability

Diagnostic-grade quality

Production according to diagnostic requirements

- ISO 13485 and ISO 9001 certifications
- Exceptionally pure enzymes and antibodies

OEM and private label

Products tailor-made to your needs

- Individual formulation, packaging and labeling
- Choice of high-quality molecular biology kits or enzymes

Lyophilization Service

Innovative lyophilization procedure

- Freeze-dried PCR enzymes and mixes
- · Custom lyophilization and LYO-ready enzymes and mixes

Contract manufacturing

Large-scale enzyme and antibody production

- Production in E. coli and eukaryotic cells
- Manufacturing and Fermentation up to 100 liter
- Purification
- Characterization
- Formulation
- Quality Control
- Assay Development
- Stability Analysis
- Dispensing and packaging

Get in touch with biotechrabbit and start your project www.biotechrabbit.com



About biotechrabbit

biotechrabbit GmbH, a German pioneering biotechnology manufacturer dedicated to producing high-quality enzymes and monoclonal antibodies tailored to meet the specific needs of diagnostic companies around the globe. As a leader in protein manufacturing, we offer private label and Original Equipment Manufacturer (OEM) solutions that empower our partners to deliver precise, reliable, and innovative diagnostic products.

Who We Are

At biotechrabbit, we harness the power of advanced biotechnology to develop and produce a wide range of proteins, with a specialization towards enzymes and antibodies for various molecular applications.

Our team of experts, equipped with the latest technology and extensive industry experience, is committed to advancing scientific progress and enhancing the accuracy and efficiency of diagnostic testing.

Our Mission

Our mission is to provide superior and high quality molecular reagents to enhance the capabilities of life science and diagnostic companies, enabling them to achieve breakthroughs in molecular applications such as medical diagnostics. We strive to support our partners in delivering consistent large-scale batches and novel developments, ultimately improving patient care and environmental research.

Innovation and Research

Innovation drives us forward. We invest heavily in research and development to stay at the forefront of enzyme technology. By fostering a culture of continuous improvement and scientific exploration, we can offer cutting-edge proteins, enzymes and formulations that keep pace with the evolving needs of the diagnostic industry.









About biotechrabbit







Partner with Us

Quality and Compliance

Quality is the cornerstone of our operations. biotechrabbit adheres to stringent quality control measures and ISO 13485/9001 regulatory standards to ensure that our products meet the highest purity, activity, and stability levels. Our state-of-the-art manufacturing facilities are equipped with advanced technologies that enable us to maintain rigorous quality assurance protocols throughout production.

Customized Enzyme Solutions

- Tailor-made enzymes and formulations are designed to meet your molecular assays' unique requirements.
- Extensive portfolio of enzymes and proteins, including a comprehensive list of polymerases, reverse transcriptases, proteases and ligases.

Private Label Manufacturing

- We offer a wide range of private label services that allow you to brand our high-quality enzymes as your own.
- Flexible production scales to accommodate your growing needs, ensuring consistent supply and reliability.

OEM Services

- Collaborative OEM partnerships focused on integrating our enzymes and formulations seamlessly into your molecular reagents.
- We provide end-to-end support from development to production, ensuring that our high-quality products enhance the performance of your molecular assays.

Join the numerous diagnostic companies that trust biotechrabbit as their preferred supplier. Whether you require customized formulations, private label manufacturing, or OEM services, we are dedicated to delivering exceptional value and performance.

Contact us today to learn more about how biotechrabbit can support your diagnostic innovations and help you achieve your goals. Together, we can significantly impact global health through superior diagnostic solutions.

Lyophilization

biotechrabbit is a premier biotechnology manufacturer dedicated to delivering innovative lyophilized products and services to diagnostic and life science sectors. Our expertise in lyophilization technology, developed over the last 13 years, allows us to offer state-of-the-art lyophilized molecular formulations that significantly enhance diagnostic assays' performance, reliability, and durability.

Our Innovative Freeze-Drying Technology

At biotechrabbit, we pride ourselves on our advanced lyophilization techniques that produce freeze-dried beads and cakes with exceptional characteristics designed to meet the rigorous demands of the diagnostic industry.

FAST DISSOLVING Our lyo beads and cakes are engineered to dissolve rapidly, ensuring swift reconstitution and enabling faster diagnostic results. This feature is crucial for time-sensitive diagnostic applications where speed and efficiency are paramount.

LOW HYGROSCOPICITY Our lyophilized beads have low moisture absorption properties, significantly reducing the risk of degradation and extending the product's shelf life. They can also be used in all kinds of cartridge production environments. This ensures your diagnostic reagents remain stable and effective over longer periods, even in challenging storage conditions.

TRIM SIZE Our production methods allow for precise and consistent dosing in the production process of beads to improve the accuracy of diagnostic tests. Their small size and various shapes also facilitate easier handling and integration into various diagnostic cartridge platforms, enhancing production convenience and overall workflow efficiency.

LOW ELECTROSTATIC CHARGE Our innovative manufacturing process ensures that our lyo beads exhibit no detectable electrostatic charge. This improves handling, simplifying the integration of the beads into automated systems and diagnostic platforms, streamlining the production process and improving overall workflow efficiency.







Lyophilization





Why Choose biotechrabbit?

Superior Performance, Quality and Compliance

biotechrabbit freeze-dries reagents using raw materials manufactures by biotechrabbit or provided by our customers. Our lyophilized products are manufactured in facilities that adhere to ISO:13485/9001 industry standards and regulatory requirements. We implement rigorous quality control measures at every stage of production to ensure that our lyophilized products meet the highest standards of purity, stability, and performance.

Customized Solutions

We understand that each partner has unique requirements. Our team of experts works closely with you to develop custom lyophilized products that precisely meet your specifications. Whether you need tailored formulations or specialized packaging, we provide solutions that align with your goals and enhance the effectiveness of your molecular assays.

Advanced Research and Innovation

Innovation drives our success. We invest heavily in research and development to continually advance our lyophilization technology and expand our product offerings. By staying at the forefront of scientific progress, we ensure that our customers benefit from the latest innovations and cutting-edge solutions in the field of lyophilization.

Partner with Us

Choosing biotechrabbit as your manufacturer for lyophilized products means partnering with a company committed to excellence, innovation, and customer satisfaction. Our dedicated team is here to support you at every step, from initial consultation and product development to production and delivery.

Our advanced freeze-drying technology not only enhances the quality and reliability of your diagnostic tests, leading to improved patient outcomes, but also supports global healthcare with a focus on sustainability. By reducing waste, minimizing energy consumption, and offering a longer shelf life, our environmentally friendly products contribute to a healthier planet while delivering superior diagnostic accuracy.

Contact us under OEM@biotechrabbit.com

Custom Protein Manufacturing

Our advanced facilities and experienced team enable us to produce recombinant proteins in both E. coli and eukaryotic cells, meeting the diverse needs of our clients in the diagnostic industry. At biotechrabbit, we leverage cutting-edge and robust biotechnological methods to produce a wide range of proteins for diagnostic applications. Our comprehensive production capabilities include:

E. coli EXPRESSION SYSTEMS

Efficient and cost-effective production of recombinant proteins with high yield and purity, ideal for various diagnostic assays. We also offer activity assay development and Quality Control.

EUKARYOTIC EXPRESSION SYSTEMS

Production of complex proteins in HEK, CHO and CapT cells with post-translational modifications, ensuring functionality and relevance for sophisticated applications.



Future Plans

We are committed to continuous improvement and expansion of our capabilities. As part of our dedication to supporting the biotech industry's evolving needs, we are in the process of establishing Good Manufacturing Practice (GMP) compliant manufacturing facilities. This will enable us to offer even higher levels of quality assurance and regulatory compliance, further solidifying our position as a leader in protein production for diagnostic and pharma applications.



Antibody Services

At biotechrabbit, we specialize in providing high-quality, customized antibody services to support your research needs. Our comprehensive offerings are designed to deliver precise and efficient results, ensuring the success of your scientific projects.

Custom Monoclonal Antibody Production

- · Choice of Species: We offer the production of monoclonal antibodies in mouse or rat.
- Starting Material: Natural or synthesized proteins and peptides or small molecules.
- Immunization and Hybridoma Development: Our efficient immunization methods, coupled with hybridoma technology, ensure successful antibody generation. This process includes extensive screening and subcloning to identify and select the best hybridoma clones.
- Proprietary Technology: We specialize in generating antibodies against challenging targets, such as membrane proteins, using our proprietary immunization and screening techniques.
- Timeline: The fast immunization protocol allows screening for hybridomas 5 weeks after the start of immunization.

Large-Scale Production

- High-Capacity Manufacturing: Our facilities are capable of producing up to 1 kg of antibodies, suitable for extensive research and commercial applications.
- Purification and Quality Control: Hybridoma cells are cultivated in serum-free media and the antibodies are purified from supernatant via protein A/G affinity chromatography, ensuring the highest purity standards. We offer mycoplasma-free and endotoxin-free production compliant with ISO:13485.

Additional Services

- Extensive Screening Assays: We provide a wide range of additional screening assays to identify the best antibody for your application.
- Flexible Manufacturing: Our manufacturing services are adaptable to your project requirements, in case of development, scale, lead time and quality.

Process Overview

- · Initial Consultation: Discuss your project requirements and goals with our experts
- Immunization and Screening: Efficient immunization methods and extensive screening to develophybridomas
- Production and Purification: Large-scale antibody production and purification to meet high-capacity demands
- · Quality Control: Stringent quality control focussing on purity, specificity and consistency

Selection Guide: Molecular Biology Enzymes & Proteins

Page Number	Product Name	PCR	RPA	LAMP	Reverse Transcription	High-Fidelity PCR	Long Range PCR	Transcription	Genome Amplification	NGS	Ligation	Sample Preparation	Cloning
11	Taq DNA Polymerase	٠											
17	YourTaq™ Hot-Start DNA Polymerase	٠											
17	Hot-Start Taq DNA Polymerase	٠											
21	ApStarTaq™ Hot-Start DNA Polymerase	٠											
24	24 AnyScript [™] Reverse Transcriptase			•	•								
24	AllScript™ Reverse Transcriptase	٠	•		•								
25	RevertUp™ II Reverse Transcriptase	٠		•	٠								
25	MMuLV Reverse Transcriptase	٠			•								
43	Long Range DNA Polymerase	٠					•						
44	Pfu DNA Polymerase	٠				•				•			
45	T4 DNA Ligase Rapid	٠								•	•		•
45	ExactiFi™ Hot-Start DNA Polymerase	٠				•				•			
49	LYO-ready T6 UvsX DNA Recombinase		•										
49	LYO-ready T4 UvsY Protein		•										
50	LYO-ready Exonuclease III		•						٠				
50	LYO-ready RB69 Gene 32 Protein	٠	•			•	•		٠				
51	LYO-ready T4 Gene 32 Protein	٠	•			٠	•		٠				
51	RNase Inhibitor	٠			٠	٠	•	•					
52	2 Heat-Labile Uracil-DNA Glycosylase												
52	Proteinase K											•	
53	3 T7 RNA Polymerase							•					
53	phi29 DNA Polymerase								٠				
58	Exo SAP Kit											•	٠



Selection Guide: biotechrabbit PCR Reagents

age Number	Product Name	Standard PCR	Direct-Load	Hot-Start PCR	Multiplex PCR	Reverse Transcription	Quantitative PCR	All-in-One Mix	-ower Crosscontamination	Green Dye included	-yophilized	-yo-Ready	-ong Range PCR	High-Fidelity PCR
		0,	-	-	~	-	U	1	_	0				-
12	Your Iaq [™] PCR Master Mix	•	-											
12	Your Taq ¹ PCR Master Mix Direct-Load	-	•											
13	AmPelity™ PCR Master Mix	•												
13	Green PCR Master Mix Direct-Load		•	-						•				
10				•										
18	Your Taq' Hot-Start PCR Master Mix			•										
18	Hot-Start PCR Master Mix		-	•										
19	Your Taq' Hot-Start PCR Master Mix Direct-Load		•	•										
19	Green Hot-Start PCR Master Mix Direct-Load		•	•						•				
21	Apstar raq ¹ ^m Hot-Start PCR Master Mix	•		•										
	Multiplay Hat Start DCD Master Mix													
22	1 Stop DT DCD Master Mix			•	•									
20	CADITAL M aDCD Probe Master Mix	•				•								
29	CAPITAL GEORFIODE Master Mix													
29	CAPITAL I-Step on the Master Mix					•	•		•					
30	CAPITAL GEOGRAPHIC MISSION MISSION AND A STATE OF THE ST						•		•					
31	CAPITAL Polep gran Master Mix					•								
31	CAPITAL ™ 1-Sten gBT-PCB Green Master Mix						•			•				
32	5Star™ dPCB Probe Master Mix					•				•				
32	UniFi TM 1-Step dBT-PCB Probe Master Mix													
02						•	•	•				•		
35	I YO-ready CAPITAI ™ αPCB Probe Mix											•		
35	I YO-ready CAPITAL™ gBT-PCB Probe Master Mix											•		
36	I YO-ready CAPITAL™ gPCB Probe Mix Plus								•			•		
36	I YO-ready CAPITAI ™ 1-Step gBT-PCB Probe Mix Plus								•			•		
39	YourTag™ Hot-Start PCB Master Mix, lyophilized	•									•			
39	9 Hot-Start PCB Master Mix Ivonhilized										•			
40	0 CAPITAL [™] gPCR Probe Master Mix, Ivophilized										•			
40	CAPITAL [™] 1-Step gRT-PCR Probe Master Mix. Ivophilized										•			
41	Long Range PCR Master Mix												٠	
42	Pfu PCR Master Mix													٠

Standard PCR

biotechrabbit Standard PCR products include highly purified YourTaq[™] DNA Polymerase which is optimized for high yield of amplification of up to 3 kb DNA targets, even from low copy number. The enzyme shows excellent PCR specificity and sensitivity for a broad range of amplicons. YourTaq Polymerase is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors. PCR mixes combine YourTaq, optimized buffer and high-quality dNTPs for the most demanding PCR applications. In addition, the enzyme is suitable for amplification of GC-rich templates (up to 70 %) by using 5X PCR Enhancer.

- Diagnostic-grade quality top performance for demanding application
- Convenient master mixes make your job easier
- Efficient abundant DNA ready for

downstream applications

• Resistant — direct amplification from crude

samples

Product	Main feature	Main PCR application	Simplified setup	Flexibility	Direct gel loading
Taq DNA Polymerase, 5 U/μl	MgCl ₂ supplied in a separate vial for maximum flexibility	Standard and demanding		\checkmark	
2X YourTaq™ PCR Master Mix	Premixed PCR reagents; just add template, primers and water	Routine, high- throughput	\checkmark		
2X YourTaq™ PCR Master Mix Direct-Load	Premixed PCR reagents including green-colored buffer for direct loading onto gel after the PCR	Routine, high- throughput with direct gel loading	~		~
2X AmPelify PCR Master Mix	Premixed PCR reagents; just add template, primers and water	Routine, high- throughput	~		
2X Green PCR Master Mix Direct-Load	Optimized Green PCR Master Mix for fast setup and direct loading on the gel	Routine, high- throughput with direct gel loading	~		~



Taq DNA Polymerase, 5 U/µl

Cat. no.	Size	Package content
BR0100102	500 U	 Taq DNA Polymerase, 5 U/µl
BR0100103	2,500 U	 5X Reaction Buffer
		 50 mM MgCl₂

FEATURES

- High product yields and robustness in a wide application range
- Highest quality utilized in molecular diagnostics and research
- Exceptionally pure Taq DNA Polymerase for routine and demanding PCR applications

APPLICATIONS

- Routine and applied PCR up to 3 kb
- RT-PCR
- TA cloning

DESCRIPTION

biotechrabbit[™] *Taq* DNA Polymerase is a first-choice enzyme for all routine and molecular diagnostics PCR applications. The exceptional quality and purity of the enzyme ensures highest performance that is required by the diagnostics industry and research labs.

The polymerase is suitable for standard and fast PCR applications ensuring high product yields from various templates with targets of up to 3 kb in size.

biotechrabbit *Taq* DNA Polymerase is a thermostable, highly processive $5' \rightarrow 3'$ DNA polymerase that has $5' \rightarrow 3'$ exonuclease activity and lacks $3' \rightarrow 5'$ exonuclease (proofreading) activity. The latter allows incorporation of modified nucleotides.

The enzyme also exhibits deoxynucleotidyl transferase activity that results in the addition of extra A overhang at the 3' ends of PCR products, allowing easy cloning of PCR products into vectors with T overhangs.

QUALITY CONTROL — ensuring highest performance

Functional assay

Human genomic DNA is amplified using the *Taq* DNA Polymerase and specific primers to produce a distinct band of 750 bp.

Self-priming activity

Standard PCR is carried out without primers, using the DNA Polymerase and human genomic DNA. No products were amplified.

Exonuclease assay

Linearized lambda/HindIII fragments are incubated with the DNA Polymerase in a 50 μl reaction mixture for 4 h at 37 °C. No degradation of DNA was observed.

Endonuclease assay

lambda DNA is incubated with the DNA Polymerase in a 50 μl reaction mixture for 4 h at 37 °C. No degradation of DNA was observed.

Nick activity

Supercoiled plasmid DNA is incubated with the DNA Polymerase in a 50 μ l reaction mixture for 4 h at 37 °C. No conversion of covalently closed circular DNA to nicked DNA was detected.

E. coli genomic DNA contamination assay

A sample of the denatured DNA Polymerase is analyzed with specific primers targeting the 16S rRNA gene in qPCR for the presence of contaminating *E. coli* DNA. No *E. coli* DNA was detectable.

Unit Definition

One unit is defined as the amount of *Taq* DNA Polymerase required to catalyze the incorporation of 10 nmol of dNTP into acid-insoluble form in 30 minutes at 72 °C in the presence of the reaction buffer.

2X YourTaq[™] PCR Master Mix

Cat.no.	Size	Package content
BR0102201	200 rxns of 50 µl	 2X YourTaq PCR
BR0102202	1,000 rxns of 50 µl	Master Mix
BR0102203	4,000 rxns of 50 µl	

FEATURES

- Exceptionally pure YourTaq DNA Polymerase
- · Optimized Master Mix for increased yield of amplification
- · Resistant to PCR inhibitor carry-over
- Excellent PCR specificity and sensitivity for a broad range of amplicons

APPLICATIONS

- Routine and demanding PCR amplification up to 3 kb
- Suitable for amplification of low target copy number
- TA cloning

DESCRIPTION

biotechrabbit[™] YourTaq PCR Master Mix is optimized for high yield of amplification of up to 3 kb DNA targets, even from low copy number. The mix shows excellent PCR specificity and sensitivity for a broad range of amplicons. YourTaq PCR Master Mix is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors.

The 2X YourTaq PCR Master Mix contains pure biotechrabbit YourTaq DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added. In addition, the mix is suitable for amplification of GC-rich templates (up to 70%).

2X YourTaq[™] PCR Master Mix Direct-Load

Cat.no.	Size	Package content
BR0102301	200 rxns of 50 µl	 2X YourTaq
BR0102302	1,000 rxns of 50 µl	PCR Master Mix
BR0102303	4,000 rxns of 50 µl	Direct-Load

FEATURES

- Exceptionally pure YourTaq DNA Polymerase
- Optimized Master Mix for increased yield of
 amplification and direct loading on the gel
- · Resistant to PCR inhibitor carry-over
- Excellent PCR specificity and sensitivity for a broad range of amplicons

APPLICATIONS

- Routine and demanding PCR amplification up to 3 kb
- Suitable for amplification of low target copy number
- TA cloning

DESCRIPTION

biotechrabbit[™] YourTaq PCR Master Mix Direct-Load is optimized for high yield of amplification of up to 3 kb DNA targets, even from low copy number. YourTaq Hot-Start PCR Master Mix shows excellent PCR specificity and sensitivity for a broad range of amplicons. The mix is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors.

The 2X YourTaq PCR Master Mix Direct-Load contains pure biotechrabbit YourTaq DNA Polymerase, extremely high-quality dNTPs, two dyes (blue and yellow) that separate during electrophoresis, allowing migration progress to be monitored, and sufficient buffer density for direct loading onto agarose gels. In addition, the mix is suitable for amplification of GC-rich templates (up to 70 %).



2X AmPelify PCR Master Mix

Cat.no.	Size	Package content
BR0101001	100 rxns of 50 µl	 2X AmPelify PCR
BR0101002	500 rxns of 50 µl	Master Mix
BR0101004	2,000 rxns of 50 µl	

FEATURES

- Optimized PCR Master Mix for minimal hands-on and fast setup
- Exceptionally pure *Taq* DNA Polymerase and highest quality dNTPs
- High product yield and robustness in a wide application range

APPLICATIONS

- Routine and high-throughput PCR up to 3 kb
- TA cloning

DESCRIPTION

biotechrabbit[™] 2X AmPelify PCR Master Mix is a perfect choice for a fast reaction setup that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. It is designed for routine, high-throughput PCR amplification of up to 3 kb DNA targets. The 2X AmPelify PCR Master Mix contains highly purified recombinant biotechrabbit *Taq* DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water need to be added.

biotechrabbit *Taq* DNA polymerase is a thermostable, highly processive $5' \rightarrow 3'$ DNA polymerase that has $5' \rightarrow 3'$ exonuclease activity and lacks $3' \rightarrow 5'$ exonuclease (proofreading) activity. The latter allows incorporation of modified nucleotides.

The enzyme also exhibits deoxynucleotidyltransferase activity that results in the addition of extra A overhang at the 3' ends of PCR products, allowing easy cloning of PCR products into vectors with T overhangs.

2X Green PCR Master Mix Direct-Load

Cat.no	Size	Package content
BR0100401	500 rxns of 50 µl	 2X Green PCR
BR0100402	2,000 rxns of 50 µl	Master Mix Direct-Load

FEATURES

- Optimized Green PCR Master Mix for fast setup and direct loading on the gel
- Excptionally pure Taq DNA Polymerase and highest quality dNTPs
- High product yields and robustness in a wide application range

APPLICATIONS

- · High-throughput PCR and immediate gel analysis
- Routine PCR up to 3 kb
- TA cloning

DESCRIPTION

biotechrabbit[™] Green PCR Master Mix is a perfect choice for a fast reaction setup that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. Additionally the special formulation allows reactions onto be loaded onto an agarose gel directly after amplification without a separate step for adding loading dye.

Green PCR Master Mix contains two dyes (blue and yellow) that separate during electrophoresis, allowing the migration progress to be monitored. Reactions with the Green PCR Master Mix have sufficient density for direct loading onto agarose gels. Green Reaction Buffer also allows mixtures containing the enzyme to be identified.

The Green PCR Master Mix contains highly purified recombinant biotechrabbit Taq DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template DNA, PCR primers and PCR-grade water need to be added.

NOTES



Hot-Start PCR (Antibody-mediated)

biotechrabbit Hot-start PCR products include highly purified YourTaq[™] Hot-Start DNA Polymerase which is optimized for high yield of amplification of up to 5 kb DNA targets, even from low copy number. The enzyme shows excellent PCR specificity and sensitivity for a broad range of amplicons. YourTaq Polymerase is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors. PCR mixes combine YourTaq, optimized buffer and high-quality dNTPs for the most demanding PCR applications. In addition, the enzyme is suitable for amplification of GC-rich templates (up to 70 %) by using 5X PCR Enhancer.

- Convenient master mixes including a direct-load version make your job easier
- Specific and sensitive excellent performance for a broad range of amplicons
- Efficient abundant DNA even from low copy targets
- Resistant direct amplification from crude

samples

Product	Main feature	Main PCR application	Simplified setup	Flexibility	Direct gel loading
UPstart™ <i>Taq</i> Antibody,1mg/ml	Hot-start monoclonal antibody for <i>Taq</i> DNA polymerases	High sensitivity, high specificity		\checkmark	
UPstart [™] Plus <i>Taq</i> Antibody, 1 mg/ml	Offers reliable inhibition of <i>Taq</i> DNA polymerases, ensuring precise thermal activation	High sensitivity, high specificity		~	
YourTaq [™] Hot-Start DNA Polymerase, 5 U/µl	Hot-start polymerase for complex PCR with high-GC content and resistance to PCR inhibitors and carry-over	High sensitivity, high specificity, complex PCR, colony PCR		~	
Hot-Start <i>Taq</i> DNA Polymerase, 5 U/µl	Antibody-based hot-start polymerase for improved PCR specificity and yield	High sensitivity, high specificity		~	
2X YourTaq™ Hot-Start PCR Master Mix	Antibody-based hot-start for complex PCR or high-GC content and resistance to PCR inhibitor carry-over	High sensitivity and specificity, complex PCR, colony PCR, direct gel loading	~		
2X Hot-Start PCR Master Mix	Antibody-based hot-start, optimized premixed PCR reagents	High-throughput, high specificity & sensitivity	~		
2X YourTaq [™] Hot-Start PCR Master Mix Direct-Load	Optimized Master Mix for increased yield of amplification and direct loading on the gel	High-throughput, high specificity, direct gel loading	 ✓ 		~
2X Green Hot-Start PCR Master Mix Direct-Load	Hot-Start PCR master mix with high product yield and robustness, enhanced with green buffer for direct loading onto gel after PCR	High-throughput, high specificity, direct gel loading	~		~

UPstart[™] Taq Antibody, 1 mg/ml

Cat.no.	Size	Package content
BR1200103	1 mg	 UPstart Taq Antibody, 1 mg/ml

FEATURES

- Inhibition of > 99 % Taq activity at 60 °C
- 200 ng UPstart Taq Antibody are blocking 1 U Taq DNA polymerase
- · Exceptionally pure made in cell culture
- Animal-free

APPLICATIONS

- Thermolabile inhibition of Taq DNA polymerases
- Convenient hot-start PCR setup at room temperature
- Fast polymerase activation with the first PCR denaturation step

DESCRIPTION

biotechrabbit[™] UPstart *Taq* Antibody is an ultra-pure monoclonal anti-*Taq* antibody against the active centre of *Taq* DNA polymerase. It is produced in cell culture to ensure highest quality.

The antibody can be used with highly efficient *Taq* DNA polymerases, provides an excellent method for "Hot-Start" PCR and enhances PCR specificity and sensitivity. PCR hot-start prevents the formation of primer dimers and nonspecific amplification and allows convenient PCR setup at room temperature.

In the first denaturation step of the thermal cycling, the UPstart *Taq* Antibody becomes nonfunctional and the active *Taq* DNA polymerase is released. This antibody-mediated hot-start method is significantly more convenient to use than other hot-start methods. Polymerase reactivation using this antibody is faster than with methods using chemically inhibited polymerases.

UPstart[™] Plus *Taq* Antibody, 1 mg/ml

Cat.no.	Size	Package content
BR1200203	1 mg	 UPstart Plus Taq Antibody, 1 mg/ml

FEATURES

- Inhibition of > 99 % Taq activity at 60 °C
- 200 ng UPstart *Taq* Antibody are blocking 1 U *Taq* DNA polymerase
- Exceptionally pure through recombinant production
- Animal-free

APPLICATIONS

- Offers reliable inhibition of *Taq* DNA polymerases, ensuring precise thermal activation.
- · Allows for easy PCR setup at room temperature
- Facilitates rapid activation of the polymerase at the first PCR denaturation step.

DESCRIPTION

biotechrabbit[™] UPstart[™] Plus *Taq* Antibody, 1 mg/ml is a state-of-the-art recombinant monoclonal anti-*Taq* antibody designed specifically to target the active centre of *Taq* DNA polymerase. Produced through advanced recombinant techniques, this antibody guarantees unmatched consistency and purity.

The antibody can be employed with highly efficient *Taq* DNA polymerases for superior "Hot-Start" PCR performance, enhancing both specificity and sensitivity. It prevents primer dimers and nonspecific amplification and simplifies PCR setup at room temperature.

During the initial denaturation step of PCR, the UP start[™] Plus *Taq* Antibody becomes inactive, releasing the functional *Taq* DNA polymerase. This recombinant antibodymediated hot-start method offers a more convenient and reliable alternative to other hot-start techniques, enabling quicker polymerase reactivation compared to methods using chemically modified polymerases.



YourTaq™ Hot-Start DNA Polymerase, 5 U/µl

Cat.no.	Size	Package content
BR0202101	250 U	 YourTaq Hot-Start
BR0202102	1,250 U	DNA Polymerase, 5 U/µl
BR0202103	5,000 U	 5X PCR Reaction Buffer
		 50 mM MgCl₂

FEATURES

- Exceptionally pure YourTaq Hot-Start DNA Polymerase
- Optimized buffer composition for increased yield of amplification
- · Resistant to PCR inhibitor carry-over
- Excellent PCR specificity and sensitivity for a broad range of amplicons

APPLICATIONS

- Routine and demanding PCR amplification up to 5 kb
- Suitable for amplification of low target copy number
- TA cloning

DESCRIPTION

biotechrabbit[™] YourTaq Hot-Start DNA Polymerase is optimized for high yield of amplification of up to 5 kb DNA targets, even from low copy number. The enzyme shows excellent PCR specificity and sensitivity for a broad range of amplicons. YourTaq Hot-Start DNA Polymerase is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors.

YourTaq Hot-Start DNA Polymerase is a highly pure enzyme. Together with the optimized buffer and highquality dNTPs (BR0600202) a mix is achieved for the most demanding PCR applications. In addition, the enzyme is suitable for amplification of GC-rich templates (up to 70%) pairing with 5X PCR Enhancer.

Hot-Start Taq DNA Polymerase, 5 U/µl

Cat.no	Size	Package content
BR0200102	500 U	 Hot-Start Taq DNA
BR0200103	2,500 U	Polymerase, 5 U/µl
		 5X Reaction Buffer
		 50 mM MgCl₂

FEATURES

- High PCR specificity and sensitivity
- Exceptionally pure Hot-Start Taq DNA Polymerase for sensitive PCR applications and high yields
- · Antibody-based Hot-Start for fast polymerase activation

APPLICATIONS

- Hot-start PCR up to 5 kb
- Amplification of low-copy-number targets
- RT-PCR and TA cloning

DESCRIPTION

biotechrabbit[™] Hot-Start *Taq* DNA Polymerase is a firstchoice hot-start PCR enzyme for all demanding PCR applications. The enzyme ensures high product yields with low background and without primer dimer formation and nonspecific priming.

The Hot-Start *Taq* DNA Polymerase is inactive during reaction setup due to the bound antibody which is quickly released at elevated temperatures, ensuring the enzyme is active only during PCR. There is no need for prolonged heating or denaturation steps.

The optional use of 5X PCR Enhancer improves PCR results in many cases, including impure template or low template abundance.

2X YourTaq[™] Hot-Start PCR Master Mix

Cat.no.	Size	Package content
BR0202201	200 rxns of 50 µl	 2X YourTaq Hot-Start
BR0202202	1,000 rxns of 50 µl	PCR Master Mix
BR0202203	4,000 rxns of 50 µl	

FEATURES

- Exceptionally pure YourTaq Hot-Start DNA Polymerase
- · Optimized Master Mix for increased yield of amplification
- · Resistant to PCR inhibitor carry-over
- Excellent PCR specificity and sensitivity for a broad range of amplicons

APPLICATIONS

- · Routine and demanding PCR amplification up to 5 kb
- Suitable for amplification of low target copy number
- TA cloning

DESCRIPTION

biotechrabbit[™] YourTaq[™] Hot-Start PCR Master Mix is optimized for high yield of amplification of up to 5 kb DNA targets, even from low copy number. The mix shows kexcellent PCR specificity and sensitivity for a broad range of amplicons. YourTaq Hot-Start PCR Master Mix is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors.

The 2X YourTaq Hot-Start PCR Master Mix contains pure biotechrabbit YourTaq Hot-Start DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added. In addition, the mix is suitable for amplification of GC-rich templates (up to 70 %) pairing with 5X PCR Enhancer.

2X Hot-Start PCR Master Mix

Cat.no	Size	Package content
BR0200205	200 rxns of 50 µl	 2X Hot-Start PCR
BR0200206	1,000 rxns of 50 µl	Master Mix
BR0200207	4,000 rxns of 50 µl	

FEATURES

- · Highest PCR sensitivity without prolonged reactivation
- Optimized PCR Master Mix for minimal hands-on and fast setup
- Exceptionally pure Hot-Start *Taq* DNA Polymerase and highest quality dNTPs

APPLICATIONS

- High-specificity and high-throughput hot-start PCR up to 5 kb
- Amplification of low-copy-number targets
- TA cloning

DESCRIPTION

biotechrabbit[™] Hot-Start PCR Master Mix is a perfect choice for a fast reaction setup that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. It is designed for lowbackground, high-throughput PCR of up to 5 kb DNA targets.

The 2X Hot-Start PCR Master Mix contains pure biotechrabbit Hot-Start *Taq* DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added.

The Hot-Start *Taq* DNA Polymerase is inactive during reaction setup due to the bound antibody, which is quickly released at elevated temperatures, ensuring the enzyme is active only during PCR. There is no need for prolonged heating or denaturation steps. The Hot-Start minimizes primer dimers and mispriming.

The optional use of 5X PCR Enhancer improves PCR results in many cases, including impure template or low template abundance.



2X YourTaq[™] Hot-Start PCR Master Mix Direct-Load

Cat.no.	Size	Package content
BR0202301	200 rxns of 50 µl	 2X YourTaq
BR0202302	1,000 rxns of 50 µl	Hot-Start PCR
BR0202303	4,000 rxns of 50 µl	Master Mix
		Direct-Load

FEATURES

- Exceptionally pure YourTaq Hot-Start DNA Polymerase
- Optimized Master Mix for increased yield of amplification and direct loading on the gel
- · Resistant to PCR inhibitor carry-over
- Excellent PCR specificity and sensitivity for a broad range of amplicons

APPLICATIONS

- Routine and demanding PCR amplification up to 5 kb
- · Suitable for amplification of low target copy number
- TA cloning

DESCRIPTION

biotechrabbit[™] YourTaq[™] Hot-Start PCR Master Mix Direct-Load is optimized for high yield of amplification of up to 5 kb DNA targets, even from low copy number. YourTaq Hot-Start PCR Master Mix shows excellent PCR specificity and sensitivity for a broad range of amplicons. The mix is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors.

The 2X YourTaq Hot-Start PCR Master Mix Direct-Load contains pure biotechrabbit YourTaq Hot-Start DNA Polymerase, extremely high-quality dNTPs, two dyes (blue and yellow) that separate during electrophoresis, allowing migration progress to be monitored, and sufficient buffer density for direct loading onto agarose gels. In addition, the mix is suitable for amplification of GC-rich templates (up to 70%) pairing with 5X PCR Enhancer.

2X Green Hot-Start PCR Master Mix Direct-Load

Cat.no	Size	Package content
BR0200502	500 rxns of 50 µl	 2X Green Hot-
		Start PCR Master
BR0200504	2,000 rxns of 50 µl	Mix Direct-Load

FEATURES

- Optimized Green Hot-Start PCR Master Mix for fast setup and direct loading on the gel
- Exceptionally pure Hot-Start Taq DNA Polymerase and highest quality dNTPs
- High product yields and robustness in a wide application range

APPLICATIONS

- High-throughput hot-start PCR and immediate gel analysis
- High-specificity hot-start PCR up to 5 kb
- Amplification of low-copy-number targets
- TA cloning

DESCRIPTION

biotechrabbit[™] 2X Green Hot-Start PCR Master Mix Direct-Load is a perfect choice for a fast reaction setup that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. It is designed for low-background, high-throughput PCR amplification of up to 5 kb DNA targets. Additionally, the special formulation allows reactions to be loaded directly onto gels after amplification without adding additional loading dye.

2X Green Hot-Start PCR Master Mix Direct-Load contains two dyes (blue and yellow) that separate during electrophoresis, allowing migration progress to be monitored. Reactions with 2X Green Hot-Start PCR Master Mix Direct-Load have sufficient density for direct loading onto agarose gels.

Hot-Start PCR (Aptamer-mediated)

biotechrabbit's ApStarTaq[™] Hot-Start DNA Polymerase is a highly purified enzyme, optimized for the high-yield amplification of up to 5 kb DNA targets, even from low copy numbers. ApStarTaq exhibits exceptional PCR specificity and sensitivity across a broad range of amplicons. Engineered to resist common PCR inhibitors such as blood (up to 20 %), ethanol, and humic acid, ApStarTaq ensures reliable PCR amplification from challenging DNA templates. Additionally, ApStarTaq is ideal for amplifying GC-rich templates (up to 70 %) when used with our 5X PCR Enhancer, ensuring robust and consistent performance.

Our ApStarTaq[™] Hot-Start PCR Master Mix combines this superior enzyme with an optimized buffer and high-quality dNTPs, designed to meet the demands of the most rigorous PCR applications. This mix provides a convenient, all-in-one solution for achieving high-specificity and high-sensitivity amplifications, making it an excellent choice for a variety of PCR applications.

- Convenient master mixes including a direct-load version make your job easier
- Specific and sensitive excellent performance for a broad range of amplicons
- Efficient abundant DNA even from low copy targets
- Resistant direct amplification from crude samples

Product	Main feature	Main PCR application	Simplified setup	Flexibility	Direct gel loading
ApStarTaq [™] Hot-Start DNA Polymerase, 5 U/µl	Aptamer-based hot-start for fast PCR, high specificity and yield	Fast PCR, high sensitivity and high specificity		\checkmark	
ApStarTaq [™] Hot-Start PCR Master Mix	Aptamer-based hot-start PCR, optimized premixed PCR reagents	Fast PCR, high sensitivity and high specificity	~	\checkmark	



ApStarTaq[™] Hot-Start DNA Polymerase, 5 U/µl

Cat.no.	Size	Package content
BR0201501	250 U	 ApStarTaq Hot-Start
BR0201502	1,250 U	DNA Polymerase, 5 U/µl
BR0201503	5,000 U	 5X Reaction Buffer
		 50 mM MgCl₂

FEATURES

- Aptamer-based hot-start functionality
- No activation steps needed
- High PCR specificity and sensitivity
- Exceptionally pure *Taq* DNA Polymerase for sensitive PCR applications and high yields

APPLICATIONS

- · Fast PCR reactions in endpoint and real-time analysis
- Hot-start PCR up to 5 kb
- Amplification of low-copy-number targets
- RT-PCR and TA cloning

DESCRIPTION

biotechrabbit[™] ApStarTaq Hot-Start DNA Polymerase is an aptamer based hot-start enzyme and the first-choice for fast PCR reactions. The exceptional quality and purity of the enzyme ensures highest performance, suitable for standard and fast PCR cycling in both, endpoint and real-time assays. It ensures high product yields with low background and without primer dimer formation or nonspecific priming.

The aptamer binds to *Taq* DNA Polymerase and inhibits the enzyme activity at temperatures below 45 °C. This ensures full hot-start functionality. The enzyme is released during standard PCR cycling conditions. There is no need for separate heating or denaturation steps, allowing fast PCR reactions.

2X ApStarTaq[™] Hot-Start PCR Master Mix

Cat.no.	Size	Package content
BR0201601	200 rxns of 50 µl	 2X ApStarTaq
BR0201602	1,000 rxns of 50 µl	Hot-Start PCR
BR0201603	4,000 rxns of 50 µl	Master Mix

FEATURES

- · Aptamer-based hot-start functionality
- No activation steps needed
- · High PCR specificity and sensitivity
- Exceptionally pure Polymerase for sensitive PCR applications and high yields

APPLICATIONS

- · Fast PCR reactions in endpoint and real-time analysis
- Hot-start PCR up to 5 kb
- Amplification of low-copy-number targets
- RT-PCR and TA cloning

DESCRIPTION

biotechrabbit[™] ApStarTaq Hot-Start PCR Master Mix is a perfect choice for a fast PCR reactions. It reduces the time required for pipetting and eliminates the need for buffer optimization.

The mix is optimized for low-background, high-throughput PCR of up to 5 kb DNA targets. The 2X Master Mix contains pure biotechrabbit ApStarTaq Hot-Start DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR-grade water are added.

The aptamer binds to *Taq* DNA Polymerase and inhibits its activity at temperatures below 45 °C. This ensures full Hot-Start functionality. The enzyme is released during standard PCR cycling conditions. There is no need for separate heating or denaturation steps, allowing fast PCR reactions. The hot-start functionality minimizes primer dimers and mispriming.

Multiplex Standard PCR

Multiplex PCR enables the amplification of multiple targets in one single PCR experiment using multiple sets of primers. The ability to reduce the number of reactions needed to test a sample for different targets helps saving time and money and therefore makes multiplex systems especially useful when large sample numbers have to be processed.

Multiplex PCR is widely used in biological and medical applications such as genotyping or different DNA testing methods in research, forensic and diagnostic laboratories. Examples for Multiplex PCR applications are genotyping (deletions, mutations, SNPs), detection and typing of pathogens or GMOs, analysis of satellite DNA (microsatellites, STRs, VNTRs), forensic analysis, food analysis, quantitative and reverse transcription PCR assays for gene expression, and more.

Multiplex Mastermixes contain all components necessary for the reaction and minimize the need for optimization thus making the development and set-up of multiplex PCR assays easy and fast.

- Optimized formulation for successful and robust performance
- Universal use suited for many types of multiplex applications
- Hot-start technology for highest specificity and sensitivity

2X Multiplex Hot-Start PCR Master Mix

Cat. no.	Size	Package content
BR0200801	50 rxns of 50 µl	 2X Multiplex Hot-
BR0200802	250 rxns of 50 µl	Start PCR Master
BR0200804	1,000 rxns of 50 µl	Mix

FEATURES

- Excellent performance and robustness in multiplex PCR
- · Optimized Master Mix for minimal hands-on and fast setup
- · Hot-start for highest sensitivity and specificity
- Exceptionally pure Hot-Start *Taq* DNA Polymerase and highest quality dNTPs

APPLICATIONS

- Fast and high-throughput multiplex PCR
- Parallel detection of multiple targets in a single assay
- Gene expression analysis, diagnostic and forensic genotyping
- Amplification of 50 bp to 2 kb targets

DESCRIPTION

biotechrabbit[™] Multiplex Hot-Start PCR Master Mix is a perfect choice for endpoint multiplex PCR. The unique buffer composition is optimized for robust simultaneous amplification of 10 or more targets from 50 bp – 2 kb in a single reaction.

The Master Mix contains pure biotechrabbit Hot-Start *Taq* DNA Polymerase, extremely high-quality dNTPs and optimized Multiplex PCR buffer; thus, only template, PCR primers and PCR Grade Water are added. The enhancers included in the mix enable efficient amplification of low abundant or GC-rich templates.

Hot-Start *Taq* DNA Polymerase is inactive during reaction setup due to the bound antibody, which is quickly released at elevated temperatures, ensuring the enzyme is active only during PCR. The hot-start efficiently minimizes primer dimers and mispriming.



Reverse Transcription and RT-PCR

biotechrabbit provides a selection of proprietary Moloney-Murine Leukemia Virus (MMuLV) and non-MMuLV reverse transcriptases with and without RNase-H activity. These exceptionally high performing reverse transciptases allow successful synthesis of cDNA with a length of up to 20 kb.

The cDNA kit shows excellent performance in first-strand cDNA synthesis for RT-PCR and qPCR, even with RNA templates having a high degree of secondary structures.

One-step RT-PCR kits provide optimized reagents, including Hot-Start *Taq* DNA Polymerase, for performing both reverse-transcription and amplification reactions in one tube.

A special RNase Inhibitor prevents RNA templates from degradation by pervasive RNases.

High performance — exceptional purity for

demanding applications

- Convenient excellent efficiency at high temperatures
- Reliable RNase Inhibitor for protecting
 your RNA

Product	Main PCR application	Main feature
AnyScript™Reverse Transcriptase	cDNA synthesis, cDNA library construction, RT-PCR, qRT-PCR	Improved and genetically engineered MMuLV Reverse Transcriptase with reduced RNase H activity and increased thermostability, superb sensitivity and specificity for fast RT reaction from low template input
AllScript™ Reverse Transcriptase	cDNA synthesis, two-step RT-PCR, RACE, qRT-PCR	Proprietary non-MMuLV based reverse transcriptase engineered for increased RNA affinity and sensitive transcription of a wide range of RNA templates including such with a high degree of secondary structures, RNase H activity specific to RNA hybridized to cDNA for improved 1-Step PCR
RevertUP [™] II Reverse Transcriptase	cDNA synthesis of ≥ 19 kb, two-step RT-PCR	A proprietary MMuLV reverse transcriptase engineered for increased temperature stability up to 60 °C, no RNase H activity, efficient cDNA synthesis of \geq 19 kb cDNA
MMuLV Reverse Transcriptase	cDNA synthesis, two-step RT-PCR	Exceptionally pure classical reverse transcriptase supplied with reaction buffer for standard applications
1-Step RT-PCR Master Mix	One-step RT-PCR	A blend of efficient thermostable reverse transcriptase and proprietary Ribonuclease Inhibitor ensures high cDNA yields; unique Hot-Start <i>Taq</i> DNA Polymerase in a mix with high-quality dNTPs and PCR enhancers allows sensitive, low-background amplification
cDNA Synthesis Kit	First-strand cDNA synthesis for RT-PCR and qPCR	Excellent yields at temperatures up to 55 °C and superior performance in demanding applications, including amplification of templates with a high degree of secondary structures

AnyScript[™] Reverse Transcriptase, 200 U/µI

Cat. no	Size	Package content
BR0401001	20,000 U	 AnyScript Reverse
		Transcriptase 200 U/ul
BR0401002	100,000 U	 5X Reverse Transcriptase
		Buffer

FEATURES

- Increased thermostability
- Exceptional performance for longer templates
- · Excellent cDNA yield

APPLICATIONS

- First-strand cDNA synthesis
- cDNA library construction
- RT-PCR and qRT-PCR

DESCRIPTION

biotechrabbit[™] AnyScript Reverse Transcriptase is an improved and genetically engineered MMuLV Reverse Transcriptase with reduced RNase H activity and increased thermostability. The enzyme is ultrapure and free of RNases and Nucleases known to spoil RT reactions; this makes AnyScript a perfect choice for first strand cDNA synthesis at higher temperature that MMuLV AnyScript is active up to 55 °C with high yield and full-length cDNA product. The enzyme is tailored to be employed in both two step and one step RT-PCR and qRT-PCR with high sensitivity and specificity.

Superb sensitivity and specificity make AnyScript an ideal enzyme for fast RT reaction from low template input in diagnostics application.

AllScript[™] Reverse Transcriptase, 4 U/µl

Cat. no.	Size	Package content
BR0400601	400 U	 AllScript Reverse
		Transcriptase, 4 U/µl
BR0400602	2,000 U	 5X Reverse Transcriptase
		Buffer

FEATURES

- Highly specific and sensitive RT-PCR
- Excellent performance in transcription of complex RNA secondary structures
- · High yields of cDNA even with targets in low copy number
- RNase H activity specific to RNA hybridized to cDNA for improved One-step PCR

APPLICATIONS

- Standard reverse transcription
- Synthesis of cDNA for cloning
- RT-PCR and qRT-PCR
- Rapid amplification of cDNA ends (RACE)
- RNA analysis by primer extension

DESCRIPTION

biotechrabbit[™] AllScript Reverse Transcriptase is a proprietary RT designed for highly specific and sensitive reverse transcription. It guarantees top performance in standard reverse transcription, synthesis of ds cDNA for cloning, RT-PCR and qRT-PCR, rapid amplification of cDNA ends (RACE) or RNA analysis by primer extension. It's high affinity to RNA allows transcription of complex RNA secondary structures and targets in low copy number, leading to high yields of cDNA.

AllScript Reverse Transcriptase is a multifunctional enzyme including RNA-dependent and ssDNA-dependent DNA polymerase, as well as RNase H activity. The RNase H activity is specific to RNA hybridized to cDNA, with no effect on pure RNA template, resulting in improved performance of subsequent PCR.



RevertUP[™] II Reverse Transcriptase, 200 U/µI

Cat. no.	Size	Package content
BR0400501	10,000 U	 RevertUP II Reverse
		Transcriptase, 200 U/µl
BR0400502	50,000 U	5X Reverse Transcriptase
		Buffer

FEATURES

- Improved performance for synthesis of long cDNAs (≥ 19 kb)
- Excellent efficiency at high temperatures up to 60 °C
- High sensitivity for cDNA synthesis from few copies of template

APPLICATIONS

- · First-strand cDNA synthesis
- · Generation of labeled cDNA
- RNA analysis by primer extension
- cDNA library construction
- RT-PCR

DESCRIPTION

biotechrabbit[™] RevertUP II Reverse Transcriptase is a proprietary MMuLV reverse transcriptase engineered by point mutations resulting in increased temperature stability up to 60 °C. This guarantees top performance with templates showing a high degree of secondary structures. The enzyme has no RNase H activity and ensures efficient synthesis of ≥ 19 kb cDNA.

RevertUP II Reverse Transcriptase is a DNA polymerase which uses RNA as a substrate and exhibits no measurable proofreading $3' \rightarrow 5'$ exonuclease function. This enzyme performs cDNA synthesis by extending a DNA primer annealed to an RNA template; it can also make copies of single-stranded DNA templates.

MMuLV Reverse Transcriptase, 200 U/µl

Cat. no.	Size	Package content
BR0400201	10,000 U	 MMuLV Reverse
BR0400202	50,000 U	Transcriptase, 200 U/µl
		 10X MMuLV RT Buffer

FEATURES

- Pure reverse transcriptase for cDNA synthesis
- High yields of first-strand cDNA
- High value for a fair price

APPLICATIONS

- First-strand cDNA synthesis
- · Generation of labeled cDNA
- RNA analysis by primer extension

DESCRIPTION

biotechrabbit[™] MMuLV Reverse Transcriptase is an exceptionally pure DNA polymerase which uses RNA as a substrate and exhibits no measurable proofreading 3'→5' exonuclease function. This enzyme performs cDNA synthesis by extending a DNA primer annealed to an RNA template; it can also make copies from a single-stranded DNA templates.

The enzyme is purified from a recombinant *E. coli* strain carrying the MMuLV reverse transcriptase gene.

1-Step RT-PCR Master Mix

Cat. no.	Size	Package content
BR0400102	100 rxns of 50 µl	2X One Step
BR0400103	500 rxns of 50 µl	Master Mix
	· · · · ·	 20X RT-RI Blend

FEATURES

- Efficient thermostable Reverse Transcriptase and RNase Inhibitor providing high cDNA yields
- Unique Hot-Start *Taq* DNA Polymerase in a mix with high-quality dNTPs
- PCR enhancers allowing sensitive low background amplification

APPLICATIONS

- One-step RT-PCR
- · Virus detection
- · Amplification of GC-rich and complex templates

DESCRIPTION

biotechrabbit[™] 1-Step RT-PCR Master Mix provides an easy and efficient way to perform both reverse transcription of RNA and PCR amplification of cDNA in one step. Only RNA template, primers and PCR Grade Water are added.

The 20X RT-RI Blend, which contains a blend of an efficient thermostable reverse transcriptase and a proprietary Ribonuclease Inhibitor, ensures high yields of cDNA.

The 2X One Step Master Mix contains unique Hot-Start *Taq* DNA Polymerase, dNTPs, MgCl₂ and stabilizers in an optimized buffer and provides high PCR product yields with minimal background even when using low-abundance and difficult templates. PCR enhancers included in the mix allow efficient amplification of complex templates including GC- or AT-rich sequences.

cDNA Synthesis Kit

Cat. no.	Size	Package content
BR0400401	10 rxns	 Reverse Transcriptase, 200U/µl
	of 20 µl	• 5X Reverse Transcriptase Buffer
BR0400403	125 rxns	 dNTP Mix
	of 20 µl	 RNase Inhibitor
BR0400404	250 rxns	 Hexamer and Oligo(dT) Primers
	of 20 µl	 PCR Grade Water

FEATURES

- Highly efficient synthesis of long cDNAs (≥ 19 kb)
- Excellent yields at temperatures up to 55 °C
- High sensitivity reverse transcription from low abundance template
- Superior performance in demanding applications, including templates with a high degree of secondary structure

APPLICATIONS

- First-strand cDNA synthesis for RT-PCR and qPCR
- Gene expression profiling
- RNA labeling and primer extension
- cDNA library construction

DESCRIPTION

biotechrabbit[™] cDNA Synthesis Kit provides superior components that ensure efficient first-strand cDNA synthesis from mRNA or total RNA templates. biotechrabbit Reverse Transcriptase enables highly efficient reverse transcription with increased thermostability.

biotechrabbit RNase Inhibitor is a potent non-competitive inhibitor of RNases. The combination of highly efficient cDNA synthesis, effective RNase inhibition and pure dNTPs allows high yields of cDNAs of more than 19 kb.

For greater application flexibility, hexamer primers, allowing all RNAs in the reaction to be used as templates, and an oligo(dT) primer, for the synthesis of cDNA from only poly(A) tailed mRNA, are included.



NULS

Quantitative Multiplex Real-Time PCR

biotechrabbit[™] CAPITAL[™] qPCR Probe, Green and 1-Step qRT-PCR Master Mixes are first-choice products for high performance real-time PCR. The mixes are designed to achieve excellent results in reaction specificity, efficiency, correlation coefficient and slope.

Intercalating dye-based mixes, such as CAPITAL qPCR Green Master Mix, allow nonspecific detection of double-stranded DNA and are faster and more cost-effective than probebased mixes. Probe-based mixes, such as CAPITAL qPCR Probe Master Mix, can be used specifically or universally.

biotechrabbit CAPITAL qPCR and qRT-PCR products can be used with the fast and standard modes as well as for multiplex PCR. The mixes are available with ROX[™] passive reference.

- High performance for early Ct values and excellent specificity
- Sensitive increased limit of detection and amplification of low copy targets
- Multiplex mono and multiplex applications
 in standard and fast cycling
- Compatible use on any real-time PCR

platform

Reference dye	Company	Instrument
	Analytica Jena	qTOWERr, qTOWER 2.0, qTOWER ³
	Applied Biosystems	7500, 7500 FAST, Viia [™] 7
	Bio-Rad	iCycler [®] , MyiQ [™] , iQ [™] 5, Opticon [™] , Opticon 2,
		Chromo4 [™] , MiniOpticon [™] , CFX96 [™] , CFX384 [™]
	BJS	Xxpress [®]
	Cepheid	Smartcycler [®]
Low	Eppendorf	Mastercycler® ep REALPLEX®, Mastercycler REALPLEX 2S
concentration	Illumina	Eco™
ROX	Fluidigm	BioMark™
	QIAGEN	Rotor-Gene® 3000, 6000, Q
	Roche Applied Science	LightCycler [®] 96, Lightcycler [®] 480, Lightcycler Nano
	Agilent	Mx4000P [®] , Mx3000P [®] , Mx3005P [®]
	Takara	Cycler Dice [™]
	Techne	Quantica [®] , PrimeQ
	Analytica Jena	qTOWER, qTOWER 2.0, qTOWER ³
	Applied Biosystems	7000, 7300, 7700, 7900, 7900HT,
		7900HT FAST, StepOne [™] , StepOnePlus™
	Cepheid	Smartcycler
High	Eppendorf	Mastercycler ep REALPLEX, Mastercycler REALPLEX 2S
concentration	Illumina	Eco™
ROX	QIAGEN	Rotor-Gene 3000, 6000, Q
	Roche Applied Science	LightCycler [®] 96, Lightcycler [®] 480, Lightcycler Nano
	Takara	Cycler Dice [™]
	Techne	Quantica [®] , PrimeQ
Fluorescein	Bio-Rad	iCycler, MyiQ, iQ 5

Instrument compatibility for passive reference dyes



4X CAPITAL[™] qPCR Probe Master Mix

Cat. no.	Size	Package content			
4X CAPITAL qF	4X CAPITAL qPCR Probe Master Mix				
BR0501401	200 rxns of 20 µl	 4X CAPITAL qPCR Probe 			
BR0501402	1,000 rxns of 20 µl	Master Mix			
BR0501403	4,000 rxns of 20 μl				
4X CAPITAL qPCR Probe Master Mix LROX					
BR0501501	200 rxns of 20 µl	 4X CAPITAL qPCR Probe 			
BR0501502	1,000 rxns of 20 µl	Master Mix LROX			
BR0501503	4,000 rxns of 20 µl				
4X CAPITAL qPCR Probe Master Mix HROX					
BR0501601	200 rxns of 20 µl	 4X CAPITAL qPCR Probe 			
BR0501602	1,000 rxns of 20 µl	Master Mix HROX			
BR0501603	4,000 rxns of 20 µl				

FEATURES

- Best in-class performance for both single and multiplex detection
- · Convenient master mix with high specificity in pathogen detection
- · Highly sensitive for low-abundance DNA targets

APPLICATIONS

- Standard and fast cycling qPCR with rapid extension rate for early Ct values
- For use on a wide range of probe technologies including Tagman[®], Molecular Beacons[®] and Scorpion[®] probes

DESCRIPTION

biotechrabbit[™] CAPITAL[™] qPCR Probe Master Mix for quantifying genomic, cDNA and viral sequences provides outstanding performance in single and multiplex qPCR. The high sensitivity provided by the mix is ideal for detection of low-abundance DNA targets in various applications, such as pathogen detection.

CAPITAL qPCR Probe Master Mix uses proprietary combination of enzyme and buffer chemistry for efficient extension and early Ct in single and multiplex qPCR. To enable the use of the kit on qPCR platforms with different reference dye concentration requirements, three kit formats are available: a one-step kit containing no ROX, as well as LROX and HROX versions containing ROX in the corresponding concentrations.

4X CAPITAL[™] 1-Step qRT-PCR Probe Master Mix

Cat. no.	Size	Package content	
4X CAPITAL 1-Step qRT-PCR Probe Master Mix			
BR0502001	200 rxns of 20 µl	4X CAPITAL 1-Step qRT-PCR	
BR0502002	1,000 rxns of 20 µl	Probe Master Mix	
BR0502003	4,000 rxns of 20 µl	 20X RTase with RNase Inhibitor 	
4X CAPITAL 1-Step qRT-PCR Probe Master Mix LROX			
BR0502101	200 rxns of 20 µl	4X CAPITAL 1-Step qRT-PCR	
BR0502102	1,000 rxns of 20 µl	Probe Master Mix LROX	
BR0502103	4,000 rxns of 20 µl	 20X RTase with RNase Inhibitor 	
4X CAPITAL 1-Step qRT-PCR Probe Master Mix HROX			
BR0502201	200 rxns of 20 µl	4X CAPITAL 1-Step qRT-PCR	
BR0502202	1,000 rxns of 20 µl	Probe Master HROX	
BR0502203	4,000 rxns of 20 µl	 20X RTase with RNase Inhibitor 	

FEATURES

- Best in-class performance for both single and multiplex detection
- · Convenient master mix for detection of low-copy pathogen targets
- High specificity and sensitivity across a wide range of sample sources

APPLICATIONS

- One step qRT-PCR from mRNA, total RNA and viral RNA targets
- For use with standard and fast qPCR platforms
- Single and multiplex qRT-PCR reactions

DESCRIPTION

biotechrabbit[™] CAPITAL[™] 1-Step qRT-PCR Probe Master Mix provides outstanding performance for real-time PCR quantification of RNA templates, including mRNA, total RNA and viral RNA from a wide range of targets. The master mix ensures high specificity and sensitivity in single and multiplex detection, making it the choice for extremely low-copy-number targets in pathogen detection.

CAPITAL 1-Step qRT-PCR Probe Master Mix uses proprietary reverse transcriptase technology and buffer chemistry for efficient cDNA synthesis and QPCR in a single tube. To enable the use of the kit on qPCR platforms with different reference dye concentration requirements, three kit formats are available: a one-step kit containing no ROX, as well as LROX and HROX versions containing ROX in the corresponding concentrations.

4X CAPITAL[™] qPCR Probe Master Mix Plus

Cat. no.	Size	Package content	
4X CAPITAL qPCR Probe Master Mix Plus			
BR0503101	200 rxns of 20 µl	 4X CAPITAL qPCR 	
BR0503102	1,000 rxns of 20 µl	Probe Master Mix UDG	
BR0503103	4,000 rxns of 20 µl		

FEATURES

- UDG activity helps prevent the amplification of carryover contamination from previous reactions.
- Optimized to detect low-abundance targets, ensuring accurate quantification even in samples with minimal genetic material.
- Pre-mixed, ready-to-use solution reduces setup time and minimizes the risk of errors.

APPLICATIONS

- Ideal for detecting and quantifying pathogen and disease-associated genes.
- Accurately measures changes in gene expression levels in response to various stimuli or treatments.
- Suitable for clinical diagnostics where reliability and precision are critical.

DESCRIPTION

biotechrabbit[™] CAPITAL[™] qPCR Probe Master Mix Plus is a high-performance solution designed for quantitative PCR (qPCR) applications that require utmost precision and reliability. This ready-to-use mix combines a robust *Taq* DNA polymerase, UDG (Uracil-DNA Glycosylase), and dUTP to provide enhanced protection against carryover contamination, ensuring the accuracy and reproducibility of your qPCR results.

The UDG enzyme is incorporated to degrade uracilcontaining DNA, a common byproduct of PCR carryover, while the inclusion of dUTP ensures that any contaminating DNA from previous PCRs is not amplified. This system not only improves the reliability of qPCR assays but also simplifies the workflow by preventing false positives.

4X CAPITAL[™] 1-Step qRT-PCR Probe Master Mix Plus

Cat. no.	Size	Package content	
4X CAPITAL 1-Step qRT-PCR Probe Master Mix Plus			
BR0502801	200 rxns of 20 µl	 4X CAPITAL qRT-PCR 	
BR0502802	1,000 rxns of 20 µl	Probe Master Mix UDG	
BR0502803	4,000 rxns of 20 µl	 20X RTase with 	
		RNase Inhibitor	

FEATURES

- Prevents carryover contamination with UDG and dUTP for reliable RNA quantification
- Integrates reverse transcription and PCR amplification in one tube, minimizing sample handling
- Maintains enzyme activity and stability throughout thermal cycling

APPLICATIONS

- Critical for comparing mRNA levels under various conditions
- Effectively identifies RNA viruses in health-related screenings
- · Supports studies on RNA roles in disease and therapy

DESCRIPTION

biotechrabbit[™] CAPITAL[™] 1-Step qRT-PCR Probe Master Mix Plus is a state-of-the-art solution optimized for quantitative reverse transcription PCR (qRT-PCR). This advanced mix integrates reverse transcriptase, *Taq* DNA polymerase, Uracil-DNA Glycosylase (UDG), and dUTP in a single, ready-to-use formulation, making it ideal for accurate RNA target detection and gene expression analysis.

The inclusion of UDG and dUTP enhances the mix's capability to prevent carryover contamination from previous PCR reactions. UDG excises uracil residues incorporated into DNA during PCR, and dUTP ensures that uracil-containing DNA is used, marking any potential contaminant for degradation. This dual approach significantly reduces the likelihood of amplifying carryover, thus improving the reliability of your results.


4X CAPITAL[™] qPCR Green Master Mix

Cat. no.	Size	Package content
4X CAPITAL qF	PCR Green Master Mix	
BR0501701	200 rxns of 20 µl	 4X CAPITAL qPCR
BR0501702	1,000 rxns of 20 µl	Green Master Mix
BR0501703	4,000 rxns of 20 µl	
4X CAPITAL qF	PCR Green Master Mix	LROX
BR0501801	200 rxns of 20 µl	 4X CAPITAL qPCR
BR0501802	1,000 rxns of 20 µl	Green Master Mix LROX
BR0501803	4,000 rxns of 20 µl	
4X CAPITAL qF	PCR Green Master Mix	HROX
BR0501901	200 rxns of 20 µl	 4X CAPITAL qPCR
BR0501902	1,000 rxns of 20 µl	Green Master Mix HROX
BR0501903	4,000 rxns of 20 µl	

FEATURES

- · Best in-class performance in a wide range of applications
- Highly specific amplification and excellent signal to noise ratio
- High sensitivity in amplification of low-abundance DNA targets with a wide range of linearity

APPLICATIONS

- Standard and fast cycling qPCR with rapid extension rate for early Ct values
- Accurate and robust gene expression analysis
- · Excellent performance in copy number variation analysis

DESCRIPTION

biotechrabbit[™] CAPITAL[™] qPCR Green Master Mix allows sensitive and specific amplification with an excellent signal to noise ratio and rapid extension rates. Extremely lowcopy-number targets can be detected with high efficiency over several logs of template concentration, while primer dimer formation is efficiently minimized.

CAPITAL qPCR Green Master Mix shows accurate and robust performance in wide range of applications, including gene expression and copy number variation analysis.

To enable the use of the kit on qPCR platforms with different reference dye concentration requirements, three kit formats are available: a one-step kit containing no ROX, as well as LROX and HROX versions containing ROX in the corresponding concentrations.

4X CAPITAL[™] 1-Step qRT-PCR Green Master Mix

Cat. no.	Size	Package content
4X CAPITAL 1	-Step qRT-PCR Gree	en Master Mix
BR0502301	200 rxns of 20 µl	 4X CAPITAL 1-Step qRT-PCR
BR0502302	1,000 rxns of 20 µl	Green Master Mix
BR0502303	4,000 rxns of 20 µl	 20X RTase with RNase Inhibitor
4X CAPITAL 1	-Step qRT-PCR Gree	en Master Mix LROX
BR0502401	200 rxns of 20 µl	 4X CAPITAL 1-Step qRT-PCR
BR0502402	1,000 rxns of 20 µl	Green Master Mix LROX
BR0502403	4,000 rxns of 20 µl	 20X RTase with RNase Inhibitor
4X CAPITAL 1	-Step qRT-PCR Gree	en Master Mix HRox
BR0502501	200 rxns of 20 µl	 4X CAPITAL 1-Step qRT-PCR
BR0502502	1,000 rxns of 20 µl	Green Master Mix HRox
BR0502503	4,000 rxns of 20 µl	 20X RTase with RNase Inhibitor

FEATURES

- Convenient mix for quantification of RNA templates
- Sensitive and specific amplification with rapid extension rate for early Ct values
- · Excellent linearity across a wide range of RNA dilutions

APPLICATIONS

- One step qRT-PCR from mRNA, total RNA and viral RNA targets
- · For use with standard and fast qPCR platforms

DESCRIPTION

biotechrabbit[™] CAPITAL[™] 1-Step qRT-PCR Green Master Mix allows sensitive and specific cDNA synthesis and qPCR in a single tube for quantifying mRNA, total RNA and viral RNA sequences. Extremely low-copy-number targets can be detected with high efficiency over several logs of template concentration.

CAPITAL 1-Step qRT-PCR Green Master Mix uses proprietary reverse transcriptase technology and buffer chemistry for efficient cDNA synthesis and qPCR in a single tube. To enable the use of the kit on qPCR platforms with different reference dye concentration requirements, three kit formats are available: a one-step kit containing no ROX, as well as LROX and HROX versions containing ROX in the corresponding concentratios.

2X 5Star™ qPCR Probe Master Mix

Cat. no.	Size	Package content
BC0502501	100 rxns of 20 µl	 2X5Star™qPCR
		Probe Master Mix

FEATURES

- · Robust amplification in both single and multiplex quantitative PCR
- Suitable for amplification from crude lysates and detection of low-copy numbers
- High specificity and sensitivity across a wide range of sample sources
- Excellent performance in presence of PCR inhibitors
- · Compatible with a wide range of quantitative PCR platforms

APPLICATIONS

- · Standard and fast cycling for both single and multiplex qPCR
- Suitable for routine and high-throughput laboratories
- Compatible with a wide range of probe technologies including Taqman®, Molecular Beacons® and Scorpion® probes
- For gene expression analysis, genotyping, pathogen detection and quantification

DESCRIPTION

biotechrabbit[™] 2X 5Star[™] qPCR Probe Master Mix is an antibody-mediated hot-start real-time PCR mix optimized for amplification and quantification of viral DNA, cDNA and genomic DNA from a wide range of targets. The mix is tailored for high specificity and sensitivity in single and multiplex amplification; 5Star[™] qPCR Mix is an excellent choice for the amplification of crude lysates, templates carrying PCR inhibitors and targets with low copy numbers.

5Star[™] qPCR Probe Master Mix uses a proprietary enzyme and buffer formulation suitable for fast extension and multiplexing in challenging PCR reactions.

5X UniFi™ 1-Step qRT-PCR Probe Master Mix

Cat. no.	Size	Package content
BR0504601	250 rxns of 20 µl	∙ 5X UniFi™
		1-Step gRT-PCR
BR0504602	1,000 rxns of 20 µl	Probe Master Mix

FEATURES

- · Best in-class performance for both single and multiplex detection
- Convenient all-in-one master mix
- High specificity and sensitivity across a wide range of sample sources

APPLICATIONS

- One step qRT-PCR from mRNA, total RNA and viral RNA targets
- For use with standard and fast qPCR platforms
- · Single and multiplex qRT-PCR reactions

DESCRIPTION

biotechrabbit[™] UniFi[™] 1-Step qRT-PCR Probe Master Mix provides outstanding performance for real-time PCR quantification of RNA templates, including mRNA, total RNA and viral RNA from a wide range of targets. The master mix ensures high specificity and sensitivity in single and multiplex detection, making it the choice for extremely low-copy-number targets in pathogen detection.

The mix is provided as a 5X concentrated master mix offering flexibility in pipetting schemes for template input in various PCR applications.

UniFi 1-Step qRT-PCR Probe Master Mix uses proprietary reverse transcriptase technology and buffer chemistry for efficient cDNA synthesis and qPCR in a single tube.



NULS

LYO-ready Master Mixes

biotechrabbit[™] LYO-ready products are optimized for airdrying and lyophilization, a technique that removes water from biological samples, enzymes, or reagents, resulting in a dry and stable product format that offers exceptional longevity and functionality.

With over a decade of experience in freeze-drying, biotechrabbit[™] offers specialized LYO-ready formulations for enzymes and reagents that ensure stability is not compromised during the lyophilization process.

Utilizing biotechrabbit[™] LYO-ready products for manufacturing of stable freeze-dried reagents with a highly extended shelf life, eliminates the need for a continuous cold chain during shipping and storage. This guarantees reduced-cost while increasing the sustainability.

biotechrabbit[™] LYO-ready master mixes are available in various concentrations to provide flexibility for custom production of single and multipe reactions in formats of beads and cakes, suitable for high-throughput and point-of-care cartridges.

- Optimal Stability for Prolonged Storage
- Cost-Efficient Logistics
- Effortless Reaction Setup
- Versatility Tailored to Your Needs

Product	Main feature	Main PCR application	IVD Development	UDG Treatment
14X LYO-ready CAPITAL™ qPCR Probe Mix	LYO-ready master mix including high-quality excipients	qPCR	\checkmark	
10X LYO-ready CAPITAL™ 1-Step qRT-PCR Probe Mix	LYO-ready master mix including high-quality excipients	qRT-PCR	~	
4X LYO-ready CAPITAL™ qPCR Probe Mix Plus	LYO-ready master mix including high-quality excipients and UDG	qPCR	 ✓ 	\checkmark
4X LYO-ready CAPITAL [™] 1-Step qRT-PCR Probe Mix Plus	LYO-ready master mix including high-quality excipients and UDG	qRT-PCR	\checkmark	~



14X LYO-ready CAPITAL[™] qPCR Probe Mix

Cat. no.	Size	Package content
BC0504305	500 rxns of 20 µl	 14X LYO-ready CAPITAL
		qPCR Probe Master Mix
		 4X qPCR Probe
		Reconstitution Buffer

FEATURES

- Freeze-dryable formulation including excipients and stabilizers
- Best-in-class performance in both single and multiplex quantitative amplification
- Convenient master mix for low-copy number detection
- · High specificity and sensitivity

APPLICATIONS

- For use with standard and fast qPCR platforms
- Single and multiplex qRT-PCR amplification
- Compatible with a wide range of probe technologies including Taqman[®], Molecular Beacons[®] and Scorpion[®] probes

DESCRIPTION

biotechrabbit[™] 14X LYO-ready CAPITAL[™] qPCR Probe Master Mix is a ready-to-use and concentrated formulation of all enzymes, dNTPs, excipients, and stabilizers for producing home-made lyophilizates with the option to include primers and probes. The time-effective lyophilization is possible for high and low volumes of the master mix without the collapse of the lyophilizate. The resulting lyophilizate is highly resistant to humidity and is swiftly dissolved in the provided reconstitution buffer.

The master mix is optimized for real-time PCR quantification of viral DNA, cDNA and genomic DNA from a wide range of targets. The mix ensures high specificity and sensitivity in single and multiplex amplification, making it the best choice for extremely low-copy-number targets in pathogen detection.

CAPITAL qPCR Probe Master Mix uses a proprietary enzyme and buffer formulation suitable for fast extension and multiplexing in challenging PCR reactions.

10X LYO-ready CAPITAL[™] 1-Step qRT-PCR Probe Mix

Cat. no.	Size	Package content
10X LYO-ready	CAPITAL 1-Step qR	T-PCR Probe Mix
BR0502901	500 rxns of 20 µl	 10X LYO-ready CAPITAL
		1-Step qRT-PCR Probe
		Master Mix
BR0502903	5,000 rxns of 20 µl	 5X qRT-PCR Probe
		Reconstitution Buffer
10X LYO-ready	CAPITAL 1-Step qR	T-PCR Probe Mix
BR0503001	500 rxns of 20 µl	 10X LYO-ready CAPITAL
		1-Step qRT-PCR Probe
		Master Mix
BR0503003	5,000 rxns of 20 µl	 2X qRT-PCR Probe
		Reconstitution Buffer

FEATURES

- Freeze-dryable formulation including excipients and stabilizers
- Best-in-class performance in both single and multiplex quantitative amplification
- Convenient master mix for low-copy number detection
- High specificity and sensitivity

APPLICATIONS

- One-step amplification of mRNA, total RNA and viral RNA targets
- · For use with standard and fast qPCR platforms
- Single and multiplex qRT-PCR amplification

DESCRIPTION

biotechrabbit[™] 10X LYO-ready CAPITAL[™] 1-Step qRT-PCR Probe Mix is a ready-to-use and concentrated formulation of all enzymes, dNTPs, excipients, and stabilizers for producing home-made lyophilizates with the option to include primers and probes. The time-effective lyophilization is possible for high and low volumes of the master mix without the collapse of the lyophilizate. The resulting lyophilizate is highly resistant to humidity and is swiftly dissolved in the provided reconstitution buffer. The reconstitution buffer is available in two different concentrations.

The master mix is optimized for real-time PCR quantification of RNA templates, including mRNA, total RNA, and viral RNA from a wide range of targets. The mix ensures high specificity and sensitivity in single and multiplex amplification, making it the best choice for extremely low-copy-number targets in pathogen detection.

CAPITAL qRT-PCR Probe Mix uses proprietary formulation for efficient RNA transcription and qPCR in a single tube.

4X LYO-ready CAPITAL[™] qPCR Probe Mix Plus

Cat. no.	Size	Package content
BR0505201	100 rxns of 20 µl	 4X LYO-ready CAPITAL[™] qPCR Probe Master Mix Plus
BR0505202	500 rxns of 20 µl	 4X qPCR Probe Reconstitution Buffer

FEATURES

- Freeze-dryable formulation including excipients and stabilizers
- Convenient master mix
- · Minimized risk of carryover contamination

APPLICATIONS

- · PCR amplification of viral DNA, cDNA and genomic DNA targets
- For use with standard and fast qPCR platforms
- · Single and multiplex qPCR amplification

DESCRIPTION

biotechrabbit[™] 4X LYO-ready CAPITAL[™] qPCR Probe Master Mix Plus, is a ready-to-use four times concentrated formulation including the adjusted concentration of heatlabile Uracil DNA Glycosylase (HL-UDG) minimizing the risk of carryover contamination. The mix includes all enzymes, dNTPs, excipients, and stabilizers for producing homemade lyophilizates with the option to include primers and probes. The time-effective lyophilization is possible for high and low volumes of the master mix without the collapse of the lyophilizate. The resulting lyophilizate is highly resistant to humidity and is swiftly dissolved in the provided reconstitution buffer.

The master mix is optimized for real-time PCR quantification of viral DNA, cDNA and genomic DNA from a wide range of targets. The mix ensures high specificity and sensitivity in single and multiplex amplification, making it the best choice for extremely low-copy-number targets in pathogen detection.

4X LYO-ready CAPITAL qPCR Probe Mix Plus uses proprietary enzyme and buffer formulation suitable for fast extension and multiplexing in challenging PCR reactions.

4X LYO-ready CAPITAL[™] 1-Step qRT-PCR Probe Mix Plus

Cat. no.	Size	Package content
BR0505301	100 rxns of 20 µl	 4X LYO-ready CAPITAL[™] 1-Step qRT-PCR Probe Master
BR0505302	500 rxns of 20 μl	Mix Plus 4X qRT-PCR Probe Reconstitution Buffer

FEATURES

- · Freeze-dryable formulation including excipients and stabilizers
- Convenient master mix
- Minimized risk of carryover contamination

APPLICATIONS

- One-step qRT-PCR amplification of mRNA, total RNA and viral RNA targets
- · For use with standard and fast qPCR platforms
- Single and multiplex qPCR amplification

DESCRIPTION

biotechrabbit[™] 4X LYO-ready CAPITAL[™] 1-Step qRT-PCR Probe Master Mix Plus is a ready-to-use four times concentrated formulation including the adjusted concentration of heat-labile Uracil DNA Glycosylase (HL-UDG) minimizing the risk of carryover contamination. The mix includes all enzymes, dNTPs, excipients, and stabilizers for producing home-made lyophilizates with the option to include primers and probes.

The time-effective lyophilization is possible for high and low volumes of the master mix without the collapse of the lyophilizate. The resulting lyophilizate is highly resistant to humidity and is swiftly dissolved in the provided reconstitution buffer. The master mix is optimized for real-time PCR quantification of RNA templates, including mRNA, total RNA, and viral RNA from a wide range of targets. The mix ensures high specificity and sensitivity in single and multiplex amplification, making it the best choice for extremely low-copy-number targets in pathogen detection.

4X LYO-ready CAPITAL 1-Step qRT-PCR Probe Mix Plus uses proprietary formulation for efficient RNA transcription and qPCR in a single tube.



NULS

Lyophilized Polymerases and Master Mixes

biotechrabbit[™] Lyo products combine an innovativ lyophilization process with high-quality enzymes and well established buffer or master mix compositions. Top performance in molecular biology or diagnostics applications is ensured. Lyo products are proven to show the same high activity as their liquid equivalents.

The freeze-dried enzymes avoid the need for a cool chain, leading to considerable cost savings during shipment and storage.

Covering the whole workflow from enzyme production, master mix development and lyophilization, biotechrabbit is the partner of choice for supplying molecular biology enzymes, reagents and mixes to diagnostics and life sciences.

- Diagnostic-grade quality top performance for demanding applications
- Economic no need for cooling during shipment and storage
- Robust innovative formulation for highly stable lyophilizates

Hot-Start Taq DNA Polymerase, 5 U/µl, lyophilized

Cat.no	Size	Package content
BR0201202	500 U	 Lyo Hot-Start Taq DNA
		Polymerase, 5 U/µl
BR0201203	2,500 U	Tag Reconstitution Buffer

FEATURES

- High PCR specificity and sensitivity
- Exceptionally pure Hot-Start *Taq* DNA Polymerase for sensitive PCR applications and high yields
- Antibody-based Hot-Start for fast polymerase activation

APPLICATIONS

- Hot-Start PCR up to 5 kb
- Ambient shipment and room-temperature storage
- Amplification of low-copy-number targets
- RT-PCR and TA cloning

DESCRIPTION

biotechrabbit[™] Hot-Start *Taq* DNA Polymerase, lyophilized is a freeze-dried version of its well-established liquid equivalent. The stabilized format allows shipment and storage without cooling. The provided reconstitution buffer ensures swift reconstitution and optimal reaction conditions for the enzyme.

Hot-Start *Taq* DNA Polymerase is a first–choice hot–start PCR enzyme for all demanding PCR applications. The enzyme ensures high product yields with low background and without primer dimer formation and nonspecific priming.

The Hot-Start *Taq* DNA Polymerase is inactive during reaction setup due to the bound antibody which is quickly released at elevated temperatures, ensuring the enzyme is active only during PCR. There is no need for prolonged heating or denaturation steps.

The optional use of 5X PCR Enhancer improves PCR results in many cases, including impure template or low template abundance.



2X YourTaq[™] Hot-Start PCR Master Mix, lyophilized

Cat. no.	Size	Package content
BR0202401	50 rxns	 2X Lyo YourTaq Hot-Start
	of 50 µl	PCR Master Mix
BR0202403	200 rxns	PCR Mix Reconstitution Buffer
	of 50 µl	

FEATURES

- Stable enzyme and mix for ambient shipment and room temperature storage
- Exceptionally pure YourTaq Hot-Start DNA Polymerase
- Optimized Master Mix for increased yield of amplification
- · Resistant to PCR inhibitor carry-over
- Excellent PCR specificity and sensitivity for a broad range of amplicons

APPLICATIONS

- Routine and demanding PCR amplification up to 5 kb
- · Suitable for amplification of low target copy number
- TA cloning

DESCRIPTION

biotechrabbit[™] lyophilized YourTaq[™] Hot-Start PCR Master Mix is a freeze-dried version of the well-established liquid equivalent. The stabilized format allows shipment and storage without cooling. The Master Mix is optimized for high yield of amplification of up to 5 kb DNA targets, even from low copy number. YourTaq Hot-Start PCR Master Mix shows excellent PCR specificity and sensitivity for a broad range of amplicons. The mix is resistant to PCR inhibitors, such as blood (up to 20 %), Ethanol or humic acid enabling PCR amplification from DNA templates with carry-over of PCR-inhibitors.

The 2X YourTaq Hot-Start PCR Master Mix contains pure biotechrabbit YourTaq Hot-Start DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added. In addition, the mix is suitable for amplification of GC-rich templates (up to 70 %) pairing with 5X PCR Enhancer.

2X Hot-Start PCR Master Mix, lyophilized

Cat. no.	Size	Packge content
BR0201101	50 rxns	 2X Lyo Hot-Start
	of 50 µl	PCR Master Mix
BR0201103	200 rxns	 PCR Mix Reconstitution Buffer
	of 50 µl	

FEATURES

- Room temperature stable enzymes and mixes
- Exceptionally pure Taq DNA Polymerase
- Hot-Start for high PCR specificity and sensitivity
- · Optimized Master Mix for fast setup

APPLICATIONS

- · Ambient shipment and room temperature storage
- Routine and demanding PCR applications
- PCR amplification up to 5 kb
- TA cloning

DESCRIPTION

biotechrabbit[™] lyophilized Hot-Start PCR Master Mix is a freeze dried version of the well-established liquid equivalent. The stabilized format allows shipment and storage without cooling. The Master Mix is a perfect choice for a fast reaction setup that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. It is designed for low-background, high-throughput PCR of up to 5 kb DNA targets.

The 2X Hot-Start PCR Master Mix contains pure biotechrabbit Hot-Start *Taq* DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added.

The Hot-Start *Taq* DNA Polymerase is inactive during reaction setup due to the bound antibody, which is quickly released at elevated temperatures, ensuring the enzyme is active only during PCR. There is no need for prolonged heating or denaturation steps. The Hot-Start minimizes primer dimers and mispriming.

4X CAPITAL[™] qPCR Probe Master Mix, lyophilized

Cat. no.	Size	Package content
BR0502601	200 rxns	 Lyo CAPITAL qPCR Probe
	of 20 µl	Master Mix
		 4X qPCR Probe
		Reconstitution Buffer

FEATURES

- Stable enzyme and mix for ambient shipment and room temperature storage
- Best in-class performance for both single and multiplex detection
- Convenient master mix with high specificity in pathogen detection
- Highly sensitive for low-abundance DNA targets

APPLICATIONS

- Standard and fast cycling qPCR with rapid extension rate for early Ct values
- For use on a wide range of probe technologies including Taqman[®], Molecular Beacons[®] and Scorpion[®] probes

DESCRIPTION

biotechrabbit[™] lyophilized CAPITAL[™] qPCR Probe Master Mix is a freeze-dried version of the well-established liquid equivalent. The stabilized format allows shipment and storage without cooling. The master mix is optimized for quantifying genomic, cDNA and viral sequences provides outstanding performance in single and multiplex qPCR. The high sensitivity provided by the mix is ideal for detection of low-abundance DNA targets in various applications, such as pathogen detection.

CAPITAL qPCR Probe Master Mix uses proprietary combination of enzyme and buffer chemistry for efficient extension and early Ct in single and multiplex qPCR.

5X CAPITAL[™] 1-Step qRT-PCR Probe Master Mix, lyophilized

Cat. no.	Size	Package content
BC0503202	200 rxns	 Lyo CAPITAL qRT-PCR
	of 20 µl	Probe Master Mix
		 5X qRT-PCR Probe
		Reconstitution Buffer

FEATURES

- Stable enzyme and mix for ambient shipment and room temperature storage
- Best in-class performance for both single and multiplex detection
- Convenient master mix for detection of low-copy
 pathogen targets
- High specificity and sensitivity across a wide range of sample sources

APPLICATIONS

- One step qRT-PCR from mRNA, total RNA and viral RNA targets
- For use with standard and fast qPCR platforms
- · Single and multiplex qRT-PCR reactions

DESCRIPTION

biotechrabbit[™] lyophilized CAPITAL[™] qRT-PCR Probe Master Mix is a freeze-dried version of the well-established liquid equivalent. The stabilized format allows shipment and storage without cooling. The master mix is optimized for real-time PCR quantification of RNA templates, including mRNA, total RNA and viral RNA from a wide range of targets. The mix ensures high specificity and sensitivity in single and multiplex detection, making it the choice for extremely low-copy-number targets in pathogen detection.

CAPITAL qRT-PCR Probe Master Mix uses proprietary reverse transcriptase technology and buffer chemistry for efficient cDNA synthesis and qPCR in a single tube.



NULS

Long-Range PCR, High-Fidelity and NGS

Pfu DNA Polymerase provides PCR accuracy that is approximately 10 times higher than *Taq* DNA Polymerase, and is suitable for demanding cloning, sequencing and expression applications.

Long-range PCR kits provide an optimized combination of *Taq* DNA polymerase and a proofreading DNA polymerase for longer PCR products with higher accuracy than is possible with the *Taq* enzyme alone. Long-range enzymes amplify templates of up to 30–40 kb with high fidelity.

biotechrabbit's NGS product section offers a comprehensive suite of solutions meticulously designed to enhance every step of the Next-Generation Sequencing (NGS) workflow. From library preparation to amplification, our products are trusted for their reliability and performance. With our A-tailing and Blunting Kits, along with high-fidelity Pfu Polymerase and versatile T4 DNA Polymerase, we ensure optimal results in genomics and beyond. Optimized — excellent enzyme performance

for demanding applications

- Accurate high-fidelity polymerases
- Reliable polymerases for large amplicons

Product		Main PCR application	Fidelity*	Amplicon length up to	Simplified setup
Long Range DNA Polymerase, 2.5 U/µl	DNA polymerases for long-range and high-fidelity PCR	Long-range, GC-rich templates	~4X higher	40 kb	
2X Long Range PCR Master Mix	Includes DNA polymerase and reagents for long-range PCR	High-throughput, long- and GC-rich templates	~4X higher	40 kb	~
<i>Pfu</i> DNA Polymerase, 2.5 U/μl	Ten times more accurate than Taq DNA Polymerase	High-fidelity, demanding PCR	~10X higher	4 kb	
2X <i>Pfu</i> PCR Master Mix	Includes <i>Pfu</i> DNA Polymerase and high-fidelity PCR reagents	High-throughput, high-fidelity	~10X higher	4 kb	\checkmark



Long Range DNA Polymerase, 2.5 U/µl

Cat. no.	Size	Package content
BR0300301	100 U	 Long Range PCR Enzyme
		Mix, 2.5 U/μΙ
BR0300302	500 U	 10X Long Range Reaction Buffer
		 5X PCR Enhancer

FEATURES

- High-productivity, long-range PCR
- Increased fidelity for accurate amplification of GC-rich templates
- · Polymerase mix for high yield and short cycle times

APPLICATIONS

- Long-range PCR up to 40 kb
- Amplification of GC-rich templates

DESCRIPTION

biotechrabbitTM Long Range DNA Polymerase is a firstchoice for amplification of targets up to 40 kb in size with higher accuracy than *Taq* DNA polymerase.

This specially designed blend of thermophilic polymerases is well suited for amplification of targets that are GC-rich and have complex structures.

For the most demanding applications, the supplied 5X PCR Enhancer can be optionally used to improve results when using templates with GC-rich sequences and complex structures.

Long Range DNA Polymerase produces a mixture of A-tailed and blunt-end PCR products. It is advisable to blunt products before cloning into blunt-end vectors.

2X Long Range PCR Master Mix

Cat. no.	Size	Package content
BR0300401	100 rxns of 50 µl	 2X Long Range
BR0300402	500 rxns of 50 µl	PCR Master Mix
		 5X PCR Enhancer

FEATURES

- Optimized Long Range PCR Master Mix for minimal hands-on and fast setup
- Mix of pure polymerases and highest quality dNTPs for high yield and short cycle times
- Increased fidelity for accurate amplification of GC-rich templates

APPLICATIONS

- High-throughput, long-range PCR up to 40 kb
- Amplification of GC-rich templates

DESCRIPTION

biotechrabbit[™] Long Range PCR Master Mix is a perfect choice for fast reaction setup for long-range PCR that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. It is designed for amplification of targets up to 40 kb in size. The master mix works well with GC-rich templates and amplifies DNA with a higher fidelity than *Taq* DNA polymerase.

The Long Range PCR Master Mix contains a blend of thermophilic polymerases, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added.

For the most demanding applications, the supplied 5X PCR Enhancer can be optionally used to improve results when using templates with GC-rich sequences and omplex structures.

Long Range PCR Master Mix produces a mixture of A-tailed and blunt-end PCR products. It is advisable to blunt products before cloning into the blunt-end vector.

Pfu DNA Polymerase, 2.5 U/µl

Cat. no.	Size	Package content
BR0300101	100 U	 Pfu DNA Polymerase, 2.5 U/µl
BR0300102	500 U	 10X Pfu Reaction Buffer
		 5X PCR Enhancer

FEATURES

- Accurate PCR for demanding applications
- Approximately ten times higher accuracy than *Taq* DNA Polymerase
- Proof-reading for increased fidelity

APPLICATIONS

- High-fidelity PCR
- Generation of PCR products for blunt cloning
- Site directed mutagenesis

DESCRIPTION

biotechrabbitTM *Pfu* DNA Polymerase is a highly purified thermostable recombinant proofreading DNA polymerase. *Pfu* DNA Polymerase exhibits approximately 10 times higher accuracy than *Taq* DNA polymerase and amplifies targets up to 3-4 kb in size.

The enzyme catalyzes template-dependent nucletide polymerization in the 5' \rightarrow 3' direction. Additionally the 3' \rightarrow 5' exonuclease (proofreading) activity corrects nucleotide incorporation errors, thereby increasing fidelity and accuracy of DNA polymerization. The enzyme has no 5' \rightarrow 3' exonuclease activity and no detectable reverse transcriptase activity and produces blunt-end PCR products.

For the most demanding applications, the supplied 5X PCR Enhancer can be optionally used for improving results when using templates with GC-rich sequences and complex structures.

2X Pfu PCR Master Mix

Cat. no.	Size	Package content
BR0300201	100 rxns of 50 µl	 2X Pfu PCR Master Mix
BR0300202	500 rxns of 50 µl	 5X PCR Enhancer

FEATURES

- Optimized *Pfu* PCR Master Mix for minimal hands-on and fast setup
- Pure Pfu DNA Polymerase and highest quality dNTPs
- Approximately ten times higher accuracy than *Taq* DNA Polymerase for accurate PCR in demanding applications

APPLICATIONS

- High-throughput, high-fidelity PCR
- Generation of PCR products for blunt cloning

DESCRIPTION

biotechrabbit^m *Pfu* PCR Master Mix is a perfect choice for fast, high-fidelity PCR setup that reduces the time required for calculation and pipetting and eliminates the need for buffer optimization. It is designed for routine highthroughput, high-fidelity amplification of targets up to 3-4 kb in size.

The 2X *Pfu* PCR Master Mix contains *Pfu* DNA Polymerase, extremely high-quality dNTPs and optimized PCR buffer; thus, only template, PCR primers and PCR Grade Water are added. For the most demanding applications, the supplied 5X PCR Enhancer can be optionally be used to improve results when using templates with GC-rich sequences and complex structures.

Pfu DNA Polymerase exhibits approximately 10 times higher accuracy compared to *Taq* DNA polymerase. *Pfu* DNA Polymerase produces blunt-end PCR products suitable for blunt cloning.



T4 DNA Ligase Rapid, 600 U/µl

Cat. no.	Size	Package content
BR1100301	60,000 U	 T4 DNA Ligase Rapid,
	(100 rxns)	600 U/µl
BR1100302	180,000U	 5X Rapid Ligation
	(300 rxns)	Buffer

FEATURES

- Exceptionally pure high-concentration T4 DNA Ligase
- Perfect for rapid ligation

APPLICATIONS

- Blunt and sticky end DNA ligation for cloning
- · Joining of linkers or adaptors to double-stranded DNA
- · Self-circularization of linear DNA

DESCRIPTION

biotechrabbit[™] T4 DNA Ligase Rapid is an exceptionally pure, highly concentrated ligase for applications in which high enzyme concentrations are required. It is especially recommended for fast ligations.

T4 DNA Ligase Rapid is supplied with 5X Rapid Ligation Buffer containing PEG for fast 5–10 minutes ligation, ligation of low-concentration or blunt-end DNA.

T4 DNA Ligase catalyzes the formation of a phosphodiester bond between the terminal 5' phosphate and the 3' hydroxyl groups of duplex DNA or RNA. The enzyme efficiently joins blunt and cohesive ends and repairs single stranded nicks in duplex DNA, RNA or DNA–RNA hybrids.

ExactiFi[™] Hot-Start DNA Polymerase, 1 U/µl

Cat. no.	Size	Package content
BR0300501	100 U	 HiFi Hot-Start DNA Polymerase, 1U/µl
		 5X HiFi Reaction Buffer
		 dNTP Mix (10 mM each)
		 50 mM MgCl₂

FEATURES

- Exceptional fidelity, approximately 50 times higher than Taq polymerase
- Amplifies targets up to 5–10 kb in size
- Produces blunt-end PCR products
- Efficient amplification of both low- and high-complexity targets
- High resistance to inhibitors for robust performance across a variety of sample types

APPLICATIONS

- High-fidelity PCR
- Generation of PCR products for blunt cloning

DESCRIPTION

biotechrabbit^m ExactiFi^m Hot-Start DNA Polymerase is a highly purified thermostable recombinant proofreading DNA polymerase. ExactiFi exhibits approximately 50 times higher accuracy than *Taq* DNA polymerase and amplifies targets up to 5–10 kb in size.

The enzyme catalyzes template-dependent nucleotide polymerization in the $5'\rightarrow 3'$ direction. Additionally, the $3'\rightarrow 5'$ exonuclease (proofreading) activity corrects nucleotide incorporation errors, thereby increasing fidelity and accuracy of DNA polymerization. The enzyme has no $5'\rightarrow 3'$ exonuclease activity and no detectable reverse transcriptase activity, producing blunt-end PCR products.

A-Tailing Kit

Cat. no.	Size	Package content
BR1101701	100 rxns	 20X A-Tailing Mix
		 dATP (1mM)
		 10X Activation Buffer

FEATURES

- Optimized for efficient and consistent A-tailing reactions
- Includes all necessary reagents for adding a single dA to the 3' ends of DNA fragments
- · High specificity and minimal background activity
- Suitable for downstream applications like cloning, sequencing, and next-generation sequencing library preparation

APPLICATIONS

- · Preparation of DNA for cloning into T/A cloning vectors
- Suitable for sequencing applications
- · Ideal for next-generation sequencing (NGS) library preparation
- Compatible with a broad range of DNA templates

DESCRIPTION

biotechrabbit[™] A-Tailing Kit contains the essential reagents required for the addition of an adenine nucleotide to the 3' ends of DNA fragments. This kit includes a 20X A-Tailing Mix, a 1 mM dATP solution, and a 10X Activation Buffer.

The reaction is designed to be performed in a single step and ensures a highly efficient A-tailing process, with minimal risk of overhang loss. The kit components are optimized to work together, providing consistent results with high specificity.

DNA Blunting Kit

Cat. no.	Size	Package content
BR1101101	100 rxns	 20X Blunting Enzyme Mix
		 10X Blunting Buffer
		 dNTP Mix (10 mM each)

FEATURES

- Rapid blunting reaction
- Excellent performance in blunt-ended ligation
- Ready-to-use blunted product in subsequent ligation

APPLICATIONS

- · Blunting and phosphorylation of double-stranded DNA
- Preparation of restriction enzyme digested DNA or sheared DNA for blunt-ended ligation

DESCRIPTION

biotechrabbit[™] DNA Blunting Kit shows excellent performance in blunting and phosphorylation of doublestranded DNA. The kit is compatible with a wide variety of dsDNA templates such as restriction enzyme digested DNA fragments, plasmid DNA, PCR products with dA-overhang, sheared or nebulized DNA. The kit is optimized for rapid blunting reaction.



NULS

Enzymes and Proteins for Molecular Biology

biotechrabbit[™] enzymes are suitable for the most demanding molecular biology applications. Selected for specific characteristics and excellent performance, biotechrabbit enzymes are provided in exceptionally pure preparations to ensure the integrity of your reactions.

- High performance success in your
 - molecular biology applications
- Exceptional purity reliable results in

demanding reactions

Product Applications		Features		
LYO-ready T6 UvsXDNA amplification in isothermal amplificationDNA Recombinasetechniques and clonal amplification in NGS		Exceptionally pure enzyme, provided in a buffer with low glycerol content for further use in lyophilization		
LYO-ready T4 UvsYPromotes the UvsX-catalysed strand invasion and assists in the release of ssDNA binding protein complexes in isothermal amplification		Highly purified protein, stored in a low-glycerol buffer to facilitate easy lyophilization		
LYO-ready Rapid and cooperative removal of nucleotides Exonuclease III from DNA strands in isothermal amplification and DNA sequencing		Enzyme with exceptional purity for molecular diagnostics, stored in a buffer that was optimised for further use in lyophilization		
LYO-ready RB69 Gene 32 Protein	Binds ssDNA effectively to boost efficiency and fidelity of PCR or isothermal amplification	Exceltionally pure protein, provided in a buffer with low glycerol content for further use in lyphilization		
LYO-ready T4 GenessDNA binding protein for increased efficiency32 Proteinand fidelity in PCR and isothermal amplification		Protein with exceptional purity, stored in a low-glycerol buffer to facilitate easy lyophilization		
RNase Inhibitor Helps prevent RNA degradation in cDNA synthesis, <i>in vitro</i> transcription/translation reactions or RNA purification		Potent inhibitor of a wide spectrum of ribonucleases under a variety of reaction conditions used for work with RNA, highly purified and stored in an optimized buffer		
Heat Labile Uracil– DNA Glycosylase	Elimination of carry-over contamination in PCR and post-PCR analysis such as cloning and sequencing, analysis of ancient DNA	Completely and irreversibly heat inactivated without addition of agents or inhibitors, active in common PCR and RT-PCR buffers		
Proteinase K Non-specific protein degradation, nucleic acid liquid and lyophilized purification, deactivation of nucleases in enzymatic reactions		Exceptional purity suitable for the most sensitive applications		



LYO-ready T6 UvsX DNA Recombinase, 5 mg/ml

Cat. no.	Size	Package content
BR1500201	100 µg	 LYO-ready T6 UvsX DNA
BR1500202	500 µg	Recombinase
		 10X Reaction Buffer I

FEATURES

- · Low glycerol content for further use in lyophilization
- Exceptional purity suitable for the most sensitive applications
- T6UvsX DNA Recombinase excels in promoting DNA strand exchange

APPLICATIONS

- Recombinase Polymerase Amplification (RPA)
- Clonal amplification in Next Generation Sequencing (NGS)
- · Gene editing

DESCRIPTION

biotechrabbit[™] LYO-ready T6 UvsX DNA Recombinase is an enzyme originating from the T6 bacteriophage. This 44 kDa enzyme is the RecA/Rad51 ortholog of bacteriophage T6 and thus plays a crucial role in facilitating DNA recombination, a process where genetic material is exchanged or integrated between DNA molecules. Highly involved in homologous recombination, UvsX binds to single-stranded DNA regions, aiding in the accurate pairing and exchange of genetic material. This enzyme can be applied in genetic engineering, gene editing, and biotechnology, serving as a tool for manipulating DNA sequences and studying the recombination machinery.

The unique properties of this enzyme make it the best choice for employment in Recombinase Polymerase Amplification (RPA), an isothermal DNA amplification method with broad applications such as diagnostics and clonal amplification in Next Generation Sequencing (NGS). This protein is provided in a storage buffer containing 5 % glycerol, making it suitable for further use in lyophilization.

LYO-ready T4 UvsY Protein, 2 mg/ml

Cat. no.	Size	Package content
BR1500301	100 µg	 LYO-ready T4 UvsY Protein
BR1500302	500 µg	 10X Reaction Buffer I

FEATURES

- Mediator Protein for T4/T6 UvsX Recombinase
- Low glycerol content for further use in lyophilization
- · Exceptionally pure Protein

APPLICATIONS

- Recombinase Polymerase Amplification (RPA)
- Mediation and stimulation of DNA dependent ATPase activity of UvsX Recombinase
- · Promotion of strand exchange
- Facilitating the release of single stranded DNA gp32 complex

DESCRIPTION

biotechrabbit[™] LYO-ready T4 UvsY Protein plays a crucial part within the DNA recombination and the repair apparatus of the T4 bacteriophage. Originating from the T4 bacteriophage, this 16 kDa protein plays a fundamental role in expediting homologous recombination processes and enhancing the effectiveness of the DNA repair, as well as promoting the loading of the T4 pre-synaptic filament onto single-stranded DNA (ssDNA). Through its interactions with pivotal recombination proteins, such as T4 and/or T6 UvsX Recombinase, T4 UvsY emerges as an indispensable element for upholding genome integrity and enabling genetic exchange.

This protein efficiently promotes the UvsX-catalyzed strand invasion by getting involved into ssDNA binding protein (SSB) complexes. The enzyme assists the release of the SSB and uplifts the binding of UvsX. This protein is provided in a storage buffer containing 0% glycerol, making it suitable for further use in lyophilization.

LYO-ready Exonuclease III, 100 U/µl

Cat. no.	Size	Package content
BR1500101	10,000 U	 LYO-ready Exonuclease III
BR1500102	25,000 U	 10X Exo-III Reaction Buffer
BR1500103	100,000 U	

FEATURES

- Low glycerol content for further use in lyophilization
- · High exonuclease activity for efficient removal of nucleotides
- Exceptional purity suitable for the most sensitive applications

APPLICATIONS

- Recombinase Polymerase Amplification (RPA)
- Other isothermal amplification techniques (e.g., RCA)
- Cloning of PCR products
- Unidirectional nested deletions (3'→5' exonuclease activity)
- · Generation of a single-stranded template for dideoxy
- Sequencing of DNA
- Production of intermediate products for site-directed mutagenesis
- Preparation of strand-specific probes

DESCRIPTION

biotechrabbit[™] LYO-ready Exonuclease III is a highly processive exonuclease that can continuously remove multiple nucleotides without dissociating from the DNA strand. This 31 kDa enzyme removes nucleotides from the 3' end of double-stranded DNA in a stepwise manner and continues to move in the 3'→5' direction. In contrast to other exonucleases that operate sequentially at one site, Exonuclease III exhibits the ability to simultaneously target multiple DNA molecules or multiple sites within a single DNA molecule. This capability enables the enzyme to rapidly and cooperatively remove nucleotides from DNA strands.

In addition to its exonuclease activity, the enzyme also exhibits RNase H, 3' DNA phosphatase, and apurinic DNA endonuclease activities. This enzymatic activity finds practical applications in DNA sequencing, removal of unwanted DNA sequences, and the creation of precise DNA fragments for genetic engineering. biotechrabbit Exonuclease III is provided in a storage buffer containing 5% glycerol, making it suitable for further use in lyophilization.

LYO-ready RB69 Gene 32 Protein, 10 mg/ml

Cat. no.	Size	Package content
BR1500401	500 µg	 LYO-ready RB69 Gene 32 Protein
BR1500402	1,000 µg	 10X Reaction Buffer I

FEATURES

- · Low glycerol content for further use in lyophilization
- · Binds single-stranded DNA with high affinity
- Essential for enhancing molecular amplification
- · Exceptionally pure protein

APPLICATIONS

- DNA isothermal amplifications (e.g., RPA, NASBA)
- DNA Repair and Recombination
- Primer Extension and Strand Displacement
- DNA Template Stabilization
- Single-Stranded DNA Protection

DESCRIPTION

biotechrabbit[™] LYO-ready RB69 Gene 32 Protein (RB69 gp32) is a single-stranded DNA-binding protein derived from bacteriophage RB69, essential for DNA replication, repair, and recombination processes. In biotechnology, RB69 gp32 serves as a valuable tool for enhancing various applications, such as PCR and isothermal amplifications. Its ability to stabilize replication forks, facilitate DNA synthesis, and protect DNA during repair makes it useful for improving the efficiency and fidelity of DNA amplification techniques.

Additionally, RB69 gp32's involvement in recombination processes offers potential applications in genetic engineering, gene editing, and the creation of novel DNA constructs, enabling researchers to manipulate and modify genetic material with greater precision and control. This protein is provided in a storage buffer containing 5 % glycerol, making it suitable for further use in lyophilization.



LYO-ready T4 Gene 32 Protein, 10 mg/ml

Cat. no.	Size	Package content
BR1500501	500 µg	 LYO-ready T4 Gene 32 Protein
BR1500502	1,000 µg	 10X Reaction Buffer I

FEATURES

- · Binds single-stranded DNA with high affinity
- · Low glycerol content for further use in lyophilization
- Essential for DNA replication, repair, and recombination
- processes, enhancing overall amplification techniques

 Exceptionally pure protein

APPLICATIONS

- Boosts the efficiency and yield of PCR through stabilization of replication forks
- Stabilizes DNA templates
- Facilitates isothermal amplifications by safeguarding DNA synthesis during the process
- Enables precise genetic engineering, gene editing, and construction of novel DNA constructs

DESCRIPTION

biotechrabbit[™] LYO-ready T4 Gene 32 Protein is a vital single-stranded DNA-binding protein originating from bacteriophage T4, pivotal for DNA replication, repair, and recombination processes.

In biotechnology, T4 Gene 32 Protein stands out as a valuable asset, amplifying a range of applications, including PCR and isothermal amplifications. Its inherent capacity to stabilize replication forks, expedite DNA synthesis, and shield DNA during repair proves instrumental in elevating the efficiency and fidelity of DNA amplification techniques.

Moreover, T4 Gene 32 Protein's active participation in recombination processes opens avenues for applications in genetic engineering, gene editing, and the creation of innovative DNA constructs. This capability empowers researchers to wield greater precision and control over the manipulation and modification of genetic material.

Our T4 Gene 32 Protein is delivered in an optimized glycerol-free storage buffer, ensuring its suitability for further use in lyophilization, facilitating extended storage and enhancing convenience for laboratory applications.

RNase Inhibitor, 40 U/µl

Cat. no.	Size	Package content
BR0400901	2,500 U	 RNase Inhibitor, 40 U/µl
BR0400902	10,000 U	

FEATURES

- Exceptionally pure proprietary Ribonuclease Inhibitor for demanding RNA applications
- Active under variety of reaction conditions used for work
 with RNA
- Prevention of RNA from degradation by a wide range of RNases

APPLICATIONS

- In vitro transcription/translation
- cDNA synthesis
- · RNA purification and storage

DESCRIPTION

biotechrabbit[™] RNase Inhibitor is a protein that is a potent inhibitor of a wide spectrum of ribonucleases. The RNase Inhibitor helps to prevent RNA degradation in applications like cDNA synthesis, RT-PCR, *in vitro* transcription/ translation reactions or RNA purification.

The enzyme is purified from a recombinant *E. coli* strain carrying the RNase Inhibitor gene.

Heat-Labile Uracil-DNA Glycosylase, 1 U/µl

Cat. no.	Size	Package content
BR1100701	100 U	 Heat Labile Uracil-DNA
BR1100702	500 U	Glycosylase, 1 U/µl

FEATURES

- The only Uracil–DNA glycosylase that is completely and irreversibly heat inactivated
- Heat-labile without any addition of agents or inhibitors
- Active in common PCR and RT-PCR buffers

APPLICATIONS

- Eliminates carry-over contamination in PCR, RT-PCR, qPCR and RT-qPCR
- Enables downstream post-PCR analysis such as cloning and sequencing
- Analysis of ancient DNA

DESCRIPTION

biotechrabbit[™] Heat-Labile Uracil-DNA Glycosylase selectively degrades uracil-containing PCR products. After performing PCR or RT-PCR using dUTP instead of dTTP, PCR products remain intact after treatment with Heat- Labile Uracil-DNA Glycosylase, whereas contaminating DNA (i.e., not amplified) is degraded. Heat-Labile Uracil-DNA Glycosylase is completely and irreversibly inactivated by moderate heat treatment at 50 °C, allowing contamination control in RT-qPCR. The enzyme hydrolyses the N-glycosylic bond between the deoxyribose sugar and the base in uracil-containing DNA leaving an abasic (apyrimidinic) site in DNA but does not modify uracils in RNA.

Heat-Labile Uracil-DNA Glycosylase is highly active at 20–40 °C. No cofactors or divalent cations are required for activity, and the enzyme is active in most PCR and RT-PCR buffers.

Proteinase K

Cat. no.	Size	Package content		
Proteinase K				
BR1101001	1.25 ml	 Proteinase K, 		
BR1101002	10 ml	20 mg/ml		
Proteinase K, lyophilized				
BR1100901	30 mg	 Proteinase K, 		
BR1100902	150 mg	lyophilized		

FEATURES

- Exceptional purity suitable for the most sensitive applications
- PCR-grade for highest performance
- Stable lyophilizate

APPLICATIONS

- Nonspecific protein degradation
- Nucleic acid purification
- · Deactivation of nucleases in enzymatic reactions

DESCRIPTION

biotechrabbit[™] Proteinase K is an active endopeptidase that is effective with native proteins, allowing endogenous RNases and DNases to be inactivated rapidly. The robust enzyme is stable over a wide pH range (4–12.5) and remains fully active for several hours when incubated at pH 6.5–9.5. The exceptional purity of the enzyme ensures that it is ideally suited for preparing PCR templates, as it is free of RNases, DNases and DNA.

biotechrabbit Proteinase K is available as an aqueous solution or lyophilized powder.



T7 RNA Polymerase, 50 U/µl

Cat. no.	Size	Package content
BR1101201	1,000 U	 T7 RNA Polymerase
BR1101202	5,000 U	• T7 RNA Polymerase Reaction Buffer

FEATURES

- Exceptionally pure enzyme
- Rigorously tested for DNA and RNase contamination to to prevent unwanted background transcription
- · Exceptional specificity for the T7 promoter

APPLICATIONS

- in vitro RNA synthesis from DNA template
- Radiolabeled RNA probes
- Non-isotopic RNA labeling
- Preparation of RNA vaccines
- · Guide RNA for gene targeting
- mRNA for in vitro translation and micro injection
- RNA structure, processing and catalysis studies
- RNA amplification
- · Anti-sense RNA for gene expression experiment

DESCRIPTION

biotechrabbit[™] T7 RNA Polymerase is a highly specific, DNA-dependent RNA polymerase renowned for its ability to selectively transcribe RNA from DNA sequences that are under the control of T7 phage promoters.

With a molecular weight of approximately 99 kDa, this enzyme plays a crucial role in the in vitro synthesis of RNA, offering a reliable solution for researchers who require precise and efficient RNA production. The enzyme's exceptional specificity for the T7 promoter ensures that RNA transcription is not only highly regulated but also incredibly efficient, making it an indispensable tool in the generation of high-quality RNA.

The enzyme is ideally suited for a variety of applications, including the synthesis of RNA for in vitro translation, the production of labeled RNA probes, the generation of antisense RNA, and various RNA structural and functional studies. In addition to its use in basic research, T7 RNA Polymerase is also a valuable tool in synthetic biology, where it facilitates the production of RNA transcripts for advanced applications such as CRISPR guide RNA synthesis, ribozyme production, and RNA aptamer generation.

phi29 DNA Polymerase, 10 U/µl

Cat. no.	Size	Package content
BR1100101	250 U	 phi 29 DNA Polymerase, 10 U/µl
BR1100102	1,000 U	 phi29 Reaction Buffer

FEATURES

- · Huge DNA yields, accurate amplification
- Highest processivity DNA polymerase with exceptionally strong strand displacement activity
- Exceptionally pure phi29 DNA Polymerase for demanding applications

APPLICATIONS

- Rolling circle amplification (RCA)
- Multiple displacement amplification
- Whole genome amplification
- Protein-primed and RNA-primed DNA amplification

DESCRIPTION

biotechrabbit[™] phi29 DNA Polymerase is an exceptionally pure DNA polymerase for demanding applications. The enzyme is purified from a recombinant E. coli strain carrying the phi29 DNA Polymerase gene from bacteriophage phi29.

The enzyme is a highly processive DNA polymerase (up to 70,000 base insertions per binding event) with a powerful strand-displacement activity and a $3' \rightarrow 5'$ proofreading exonuclease function. phi29 DNA Polymerase proofreading activity acts preferentially on single-stranded DNA or RNA.

Therefore, to avoid primer degradation during the DNA synthesis, 3' modified (protected) primers are highly recommended.

Nucleotides

biotechrabbit[™] deoxynucleotide triphosphates are available as mixes containing all dNTPs or sets of individual dNTPs. Sets and mixes containing dUTP instead of dTTP are available for applications such as PCR carryover prevention. The outstanding purity of the nucleotides ensures excellent performance in the most demanding applications. Reproducible — exceptional purity for

excellent results

Reliable — outstanding stability for

consistent PCR results

dNTP Sets and Mixes

Cat. no.	Size	Package content	
dNTP Set, 4 × 100 mM solutions			
BR0600601	4 × 250 µl	 dATP/dCTP/dGTP/dTTP 	
BR0600602	4 × 1 ml	(100 mM each)	
dNTP Mix (25 m	M each)		
BR0600502	1 ml	 dNTP Mix (25 mM each) 	
BR0600503	5 ml		
dNTP Mix (10 mM each)			
BR0600202	1 ml	 dNTP Mix (10 mM each) 	
BR0600204	5 ml		
dNTP/dUTP Mi	x (10 mM dATP	/CTP/GTP, 30 mM dUTP)	
BR0601101	1 ml	 dNTP/dUTP Mix 	
BR0601102	5 ml		
dUTP (100 mM solution)			
BR0601201	0.75 ml	 dUTP (100 mM) 	
BR0601202	3 ml		

FEATURES

- Exceptional quality dNTPs of > 99 % purity confirmed by HPLC
- Free from DNA and PCR inhibitors
- Consistent PCR results due to outstanding dNTPs stability

APPLICATIONS

- Standard or Hot-Start PCR
- · Long-range and high-fidelity PCR
- cDNA synthesis and RT-PCR
- qPCR
- Sequencing
- DNA labeling

DESCRIPTION

biotechrabbit[™] deoxynucleotide triphosphates are firstchoice nucleotides for all PCR applications, including the most demanding, such as amplification of long targets (up to 40 kb), GC-rich templates, qPCR, cDNA synthesis, highfidelity PCR, DNA labeling and sequencing.

Advanced production technology ensures that deoxyribonucleotide triphosphates have > 99 % purity and outstanding stability, ensuring excellent performance and consistent, reliable results.

For the maximum flexibility, nucleotides are available in sets and mixes of common concentrations.



Additives and Accessories

Discover a new standard of excellence in PCR applications with biotechrabbit's premium PCR additives and accessories. Our commitment to quality and purity sets the stage for unparalleled performance in your molecular biology experiments, ensuring reproducibility, reliability, and exceptional results.

Highest-quality raw materials

• Flexibility — adjust your reaction

50 mM MgCl₂

Cat.no.	Size	Package content
BR1900101	6 ml	 50 mM MgCl₂

FEATURES

- Flexibility in PCR optimization
- Improved PCR performance

APPLICATIONS

· Optimization of PCR buffer conditions

DESCRIPTION

biotechrabbit[™] MgCl₂ is ideally suited for optimization of PCR conditions. For convenience and flexibility, it is provided as 50 mM solution.

5X PCR Enhancer

Cat.no.	Size	Package content
BR1900201	6 ml	 5X PCR Enhancer

FEATURES

- Increased specificity and efficiency
- Improved results with GC-rich templates

APPLICATIONS

- Optimize PCR performance in demanding applications
- Sensitive, low-background amplification

DESCRIPTION

biotechrabbit[™] PCR Enhancer is a unique PCR additive that improves sensitivity, resulting in increased PCR efficiency and reduced background. PCR Enhancer is ideal for difficult templates (e.g., GC-rich) and complex reactions, such as low-abundance templates and multiplex PCR.

PCR Grade Water

Cat.no.	Size	Package content
BR1900301	15 ml	PCR Grade Water
BR1900302	300 ml	

FEATURES

- Ultrapure sterile filtrated deionized water
- Prepared without the use of chemicals such as DEPC
- DNase-, RNase- and Protease-free
- Bioburden tested

APPLICATIONS

- Ready-to-use water for standard and enhanced PCR and qPCR
- Buffer preparation and enzyme dilutions
- All other molecular biology applications

DESCRIPTION

biotechrabbit[™] PCR Grade Water is ultrapure and sterile filtered deionized water prepared by a proprietary process without the use of chemicals such as DEPC (diethylpyrocarbonate). The water is free from DNase, RNase, protease and bioburdon contaminations, for use in PCR and all other molecular biology applications.

Nucleic Acid Purification

biotechrabbit[™] nucleic acid purification kits are designed to be fast and convenient to use. Purified nucleic acids are suitable for standard and demanding applications.

Universal kits provide reagents and protocols for purification from a wide range of starting materials.Mini filter based procedures allow easy handling, and proprietary buffer compositions ensure most efficient lysis, nucleic acid binding, washing and elution.

- Convenient fast and simple procedures
 to make your work easier
- Efficient proprietary buffer compositions for high yields of pure nucleic acid

Product	Applications	Starting material
DNA cleanup		
GenUP PCR/Gel Cleanup Kit	PCR product cleanup and DNA extraction from agarose gels	Amplification reactions and TAE or TBE agarose gels
GenUP Exo SAP Kit	PCR cleanup for sequencing with 100 % product recovery	Amplification reactions
Plasmid isolation		
GenUP Plasmid Kit	Universal kit for isolating high- and low-copy-number plasmids from bacterial cultures	Gram-positive and Gram-negative bacteria
Total DNA isolation		
GenUP Bacteria gDNA Kit	Universal kit for genomic DNA isolation from bacterial cultures	Gram-positive and Gram-negative bacteria
GenUP gDNA Kit	Universal kit for genomic DNA isolation from various eukaryotic starting materials	Tissues, rodent tails, paraffin-embedded samples, swabs, eukaryotic cells suspensions
GenUP Blood DNA Kit	Optimized kit for DNA isolation from blood	Fresh or frozen whole blood stabilized with EDTA or citrate
GenUP Plant DNA Kit	Optimized kit for DNA isolation from plants	Fresh, frozen or dry plant materials



Nucleic Acid Purification

- Flexible purification for a range of sample
 - types with one kit
- Reliable best results in downstream
 applications

- Optimized — no toxic β -mercaptoethanol and

no DNase for RNA preparations

• Fast - PCR cleanup in only 3 min

Product	Applications	Starting material
Total RNA isolation		
GenUP Total RNA Kit	Universal kit for total RNA isolation from various sources	Tissue, eukaryotic cell suspensions, bacteria
GenUP Micro RNA Kit	Isolation of snRNA, miRNA, siRNA, tRNA, rRNA and mRNA	Eukaryotic cells, tissue or biopsies, Gram-posi- tive and Gram-negative bacteria
GenUP Blood RNA Kit	Optimized kit for total RNA isolation from blood	Fresh or frozen whole blood stabilized with EDTA or citrate
GenUP Plant RNA Kit	Optimized kit for total RNA isolation from plants	Fresh, frozen or dry plant materials
Virus DNA and RNA isola	ation	
GenUP™ BS Virus RNA Kit	Virus RNA isolation from swabs	Swab samples
GenUP Virus DNA/RNA Kit	Optimized kit for simultaneous viral DNA and RNA isolation from eukaryotic samples	Serum, plasma and other cell-free body fluids, supernatants from cell cultures, tissues, biopsies, cell cultures, swabs, paraffin- embedded samples
Deparaffinization		
GenUP FFPE Paraffin Removal Solution	Removal of paraffin from formalin-fixed, paraffin-embedded (FFPE) sample	FFPE slices

GenUP[™] PCR/Gel Cleanup Kit

Cat. no.	Size	Package content
BR0700501	10 preps	Solutions, Mini Filters
BR0700502	50 preps	and Vials
BR0700503	250 preps	

Gel cleanup

Starting material	TAE or TBE agarose gel	
	(up to 300 mg)	
Extraction time	Approximately 20 min	
Binding capacity	> 20 µg DNA	
Typical yield	100 bp – 30 kb	
Average purity	60–90 %	

PCR cleanup

Starting material	PCR mixtures (up to 50 µl)
Extraction time	Approximately 3 min
Binding capacity	> 20 µg DNA
Typical yield	60 bp – 30 kb
Average purity	60–95 %

FEATURES

- Dual performance kit for both PCR product cleanup and DNA purification from agarose gels
- Fast and simple procedure
- High DNA recovery yields

APPLICATIONS

- · Fast purification of DNA from agarose gels
- Fast purification of products from PCR amplification reactions

DESCRIPTION

biotechrabbit[™] GenUP[™] PCR/Gel Cleanup Kit has been specially developed for a quick and easy cleanup or concentration of PCR fragments from reaction mixtures as well as extraction of DNA from both TAE and TBE agarose gels. The DNA is bound to a Mini Filter using a novel buffer, washed and then eluted in a separate tube. The purified DNA is ready to be used in all demanding molecular biology applications, including restriction digestion, ligation, sequencing, transfection into mammalian cells and *in vitro* transcription.

GenUP[™] Exo SAP Kit

Cat. no.	Size	Package content
BR0701801	100 reactions	Exol-SAP
BR0701802	500 reactions	
BR0701803	2,000 reactions	

FEATURES

- · Fast 5 min PCR cleanup without the need for spin columns
- 100 % recovery no loss of PCR product
- Simple process just add Exonuclease 1 and Shrimp Alkaline Phosphatase to PCR mix
- · Easy scalability and automation

APPLICATIONS

- PCR cleanup prior to sequencing or genotyping
- Removal of nucleotides and primers from PCR reaction

DESCRIPTION

biotechrabbit[™] GenUP[™] Exo SAP Kit is designed for quickly and easily cleaning up PCR amplification reactions, enabling their direct use in sequencing or genotyping.

The kit provides a combination of Shrimp Alkaline Phosphatase (SAP) and Exonuclease 1 (Exol) for removing nucleotides and primers. Double stranded DNA is not affected, ensuring 100 % recovery of PCR products.

Both enzymes are active in PCR buffers and completely inactivated by heating at 80 °C: no further processing is required. Shrimp Alkaline Phosphatase dephosphorylates nucleotides and primers, and the $3' \rightarrow 5'$ exonuclease activity of Exonuclease 1 degrades single-stranded primers.



GenUP™ Plasmid Kit

Cotino	Sizo	Paakaga contont	
Gal. no.	Size	Package content	
BR0700201	10 preps	Solutions, Mini Filters and	
BR0700202	50 preps	Vials	
BR0700203	250 preps		
		·	
Starting material	Starting material Bacterial cultures: 0.5–15 ml for		
high-copy-		number plasmids, low-	
copy-number plasmids or cosmids		per plasmids or cosmids	
Extraction time	Approximately 15 min		
Binding capacity	Approximately 80 µg plasmid DNA		
Typical yield	Variable: 4	Variable: 45–60 µg high-copy	
	plasmid DN	plasmid DNA is typically purified	
	from 15 ml culture, or 6–20 µg		
	from 2 ml culture		
Average purity	A ₂₆₀ /A ₂₈₀ 1.7-2.0		

FEATURES

- Universal kit for both high- and low-copy-number plasmids isolation
- Fast and simple procedure
- High yields of pure plasmid DNA suitable for all applications

APPLICATIONS

• Plasmid DNA isolation from Gram-negative and Gram-positive bacteria

DESCRIPTION

biotechrabbit[™] GenUP[™] Plasmid Kit is designed for fast and efficient plasmid purification from 0.5–15 ml bacterial suspensions in a mini format. After performing an alkaline lysis step and precipitating chromosomal DNA and bacterial proteins, plasmid DNA is bound to a Mini Filter, washed and then eluted in a low-salt buffer. Typically, yields are up to 60 µg plasmid DNA from a 15 ml bacterial suspension, depending on a plasmid copy number.

The purified plasmid DNA is ready to be used in all demanding molecular biology applications, including enzymatic reactions and sequencing.

GenUP™ Bacteria gDNA Kit

Cat. no.	Size	Package content	
BR0700701	10 preps	Solutions, Mini Filters	
BR0700702	50 preps	and Vials	
Starting material 1×10 ⁹ Gram-negative or			
	Gram-positive bacterial cells		
Extraction time	Approxima	Approximately 45 min	
Binding capacity	> 50 µg DN	> 50 µg DNA	

Typical yield Variable; approximately 35 µg

FEATURES

- Fast and simple procedure
- gDNA from up to 1 × 10⁹ bacteria
- High yields of pure DNA for demanding applications

APPLICATIONS

 Isolation of bacterial genomic DNA from Gram-positive and Gram-negative bacteria

DESCRIPTION

biotechrabbit[™] GenUP[™] Bacterial gDNA Kit has been specially developed for quick and easy purification of bacterial genomic DNA from both Gram-negative and difficult to process Gram-positive bacteria.

A combined lysozyme and proteolytic lysis steps allow efficient cell disruption. The DNA is bound to a highcapacity filter, washed and then eluted in a separate tube. The purified DNA is ready to be used in all demanding molecular biology applications, including PCR, enzymatic digestions, cloning and other.

GenUP™ gDNA Kit

		-	
Cat. no.	Size	Package content	
BR0700601	10 preps	Solutions, Mini Filters	
BR0700602	50 preps	and Vials	
BR0700603	250 preps		
Starting material Tissue samples (up to 40 mg),			
rodent tail specimens of 0.5–1 cm			
in length, eukaryotic cells			
(up to 5×10^6), formalin-fixed) ⁶), formalin-fixed	
	paraffin-embedded tissue samples		
	(FFPE, 2–4	(FFPE, 2–4 slices), buccal swabs	
Extraction time	Approximately 8 min after lysis		
Binding capacity	> 100 µg DNA		
Typical yield	Variable; approximately 65 µg		
Average purity	A ₂₆₀ /A ₂₈₀ 1.7-2.0		

FEATURES

- · Universal kit for DNA isolation from various starting materials
- Fast and simple procedure with > 100 µg gDNA filter-binding capacity
- High yields of pure DNA

APPLICATIONS

 Isolation of genomic DNA from tissues, cells, rodent tail, buccal swaps and paraffin samples

DESCRIPTION

biotechrabbit[™] GenUP[™] gDNA Kit is designed for a fast and efficient genomic DNA isolation from various sources and different amounts of starting material, such as mammalian tissues (including paraffin-embedded), buccal swabs and eukaryotic cell cultures. After a few initial procedures, the genomic DNA is bound to a Mini Filter, washed and then eluted in a low-salt buffer. High yields of pure, unshared genomic DNA are ready to be used in all demanding molecular biology applications, including enzymatic reactions and sequencing.



GenUP[™] Blood DNA Kit

Cat. no.	Size	Package content
BR0701301	10 preps	Solutions, Mini Filters
BR0701302	50 preps	and Vials
BR0701303	250 preps	
Starting material Fresh or frozen whole blood;		ozen whole blood;
stabilized with EDTA or citrate		vith EDTA or citrate
(200 μl or 400 μl)		400 μl)
Extraction time	Approximately 24 min	
Binding capacity	> 60 µg DNA	
Typical yield	Variable depending on the starting	
	material; approximately 30 µg DNA	
Average purity	A ₂₆₀ /A ₂₈₀ 1.7-2.0	

FEATURES

- · Fast and simple procedure
- Genomic gDNA from fresh and frozen, EDTA- or citrate-treated blood
- Excellent genomic DNA quality in yields of up to 30 µg

APPLICATIONS

Isolation of genomic DNA from up to 400 µl whole blood

DESCRIPTION

biotechrabbit[™] GenUP[™] Blood DNA Kit is designed for fast isolation of genomic DNA from up to 400 µl whole blood from fresh or frozen samples that have been stabilized with EDTA or citrate. After an efficient lysis step, genomic DNA is bound to a Mini Filter, washed and eluted. The isolation chemistry and extraction protocol are optimized for maximum yield. Including lysis, isolated DNA is available in approximately 24 min. The isolated DNA is suitable for a wide range of different molecular biology applications.

Protocols are available for isolating DNA from 200 μl or 400 μl whole blood samples.

The GenUP Blood DNA Kit is designed for the use with blood. For other starting material, such as cell-free body fluids (including cerebrospinal fluid, serum, plasma or urine), tissue, stool samples, buffy coat, cultured or isolated cells, swabs, dried blood spots, viruses, fungi, bacteria or parasites, please refer to the GenUP gDNA Kit (cat. no. BR0700601), GenUP Bacteria gDNA Kit (cat. no. BR0700701), GenUP Plant DNA Kit (cat. no. BR0700801) or GenUP Virus DNA/RNA Kit (cat. no. BR0701101).

GenUP[™] Plant DNA Kit

Cat. no.	Size	Package content
BR0700801	10 preps	Solutions, Mini Filters
BR0700802	50 preps	and Vials
BR0700803	250 preps	
Starting material Fresh, frozen or dried plant tissue		en or dried plant tissue
(maximur		100 mg dry weight
or 180 mg		wet weight)
Extraction time	n time Approximately 40 min	
Binding capacity	> 50 µg DNA	
Typical yield	Variable depending on the starting ma-	
	terial; approximately 5–25 µg DNA	
Average purity	A ₂₆₀ /A ₂₈₀ 1.7-2.0	

FEATURES

- Simple and efficient procedure for plant DNA isolation
- Special lysis protocols for different plant materials
- High yields of inhibitor-free DNA

APPLICATIONS

 Universal kit for isolating genomic DNA from various plant materials

DESCRIPTION

biotechrabbit[™] GenUP[™] Plant DNA Kit has been specially developed for quick and easy purification of genomic DNA from a wide variety of plant materials, including fresh, frozen or dried samples from leaves, roots, stems and flowers.

The kit includes an advanced prefiltration step to remove unlysed tissue. Subsequently, DNA is bound to a Mini Filter and is subsequently washed and eluted in a separate tube. The purified DNA is ready for use in any demanding molecular biology application, including PCR, enzymatic digestions and cloning.

This kit provides three buffers for optimized processing with different plant materials. To determine optimal lysis conditions, side-by-side preparation using the three provided protocols are prepared.

GenUP[™] Total RNA Kit

Cat. no.	Size	Package content
BR0700901	10 preps	Solutions, Mini Filters
BR0700902	50 preps	and Vials
BR0700903	250 preps	
Starting material Eukaryotic cells (5 × 10 ⁶),		
tissue samples (up to 20 mg), bacteria		
	cells (Gram-positive or Gram-negative,	
	1 × 10 ⁹)	
Extraction time	Approximately 20–40 min	
Binding capacity	100 µg RNA	
Typical yield	Yield is highly dependent on	
	sample tvp	e

FEATURES

- Fast and simple procedure
- High yields of pure RNA
- Physical removal of DNA, no DNase treatment, no toxic β-mercaptoethanol

GenUP™ Micro RNA Kit

Cat. no.	Size	Package content	
BR0701901	10 preps	Solutions, Mini Filters	
BR0701902	50 preps	and Vials	
BR0701903	250 preps		
Starting material	Eukaryotic	cells (5 ×10 ⁶)	
	Fresh or fro	ozen tissue or biopsies (up	
	to 20 mg)	to 20 mg)	
	Gram-positive and Gram-negative		
	bacteria (up to 1 × 10 ⁹)		
Extraction time	Typically 15	5–40 minutes	
Binding capacity	100 µg		
Typical yield	Depends o	Depends on the type and the amount	
	of the start	of the starting material	
Recovery rate	High rate of recovery for small RNA		
	molecules		

APPLICATIONS

 Universal kit for total RNA isolation from various sources and different amounts of starting material

DESCRIPTION

biotechrabbit[™] GenUP[™] Total RNA Kit has been specially developed for a quick and easy purification of total RNA from eukaryotic cell suspensions, tissues and biopsies, Gram-negative (e.g., *E. coli*) and Gram-positive bacteria and other sources. After few initial procedures, the RNA is bound to a filter, washed and then eluted in a separate tube. DNA is removed physically by binding to a filter without any DNAse treatment or the use of toxic β-mercaptoethanol.

The purified RNA is ready to be used in all demanding molecular biology applications, including cDNA synthesis, northern blot analysis and others.

FEATURES

- Fast and simple procedure
- Optimized binding conditions for high yields
 of small RNAs
- Physical removal of DNA, no DNase treatment, no toxic β-mercaptoethanol

APPLICATIONS

• Efficient isolation of snRNA, miRNA, siRNA, tRNA, rRNA and mRNA from various starting material

DESCRIPTION

biotechrabbit[™] GenUP[™] Micro RNA Kit is designed for high yields of small RNA molecules and total RNA, including snRNA, miRNA, siRNA, tRNA, rRNA and mRNA, without the use of highly toxic β-mercaptoethanol. After using well established filter-technology to selectively remove genomic DNA, the RNA is bound, washed and eluted from the filter membrane using RNase-free water.



GenUP[™] Blood RNA Kit

Cat. no.	Size	Package content	
BR0701401	10 preps	Solutions, Mini Filters	
BR0701402	50 preps	and Vials	
BR0701403	250 preps		
Starting material Fresh or frozen whole blood;			
stabilized with EDTA or citrate			
(0.5–1.0 ml)			
Extraction time Approximately 45 min			
Binding capacity	Binding capacity > 20 μg RNA		
Typical yield	Variable depending on the starting		
	material; 1–8 µg RNA		

FEATURES

- Fast and simple procedure
- High-quality RNA isolated from fresh and frozen, EDTA- or citrate-treated blood
- Physical removal of DNA, no DNase treatment, no toxic β-mercaptoethanol

APPLICATIONS

 Isolation of up to 8 µg total RNA from 0.5–1.0 ml whole blood

GenUP™ Plant RNA Kit

Cat. no.	Size	Package content	
BR0701501	10 preps	Solutions, Mini Filters	
BR0701502	50 preps	and Vials	
Starting material Plant material (up to 100 mg)			
Extraction time	e 30 min after homogenization		
Binding capacity	Approximately 100 µg RNA		
Typical yield	Variable depending on the starting		
	material; approximately 70 µg RNA		
Average purity	Apen/Apen 17-20		

FEATURES

- · Fast and simple procedure
- High-quality RNA isolated from a wide variety of plant samples
- Physical removal of DNA, no DNase treatment, no toxic β-mercaptoethanol

APPLICATIONS

• Isolation of total RNA from up to 100 mg plant material

DESCRIPTION

biotechrabbit[™] GenUP[™] Blood RNA Kit is designed for fast isolation of total RNA from up to 1 ml whole blood from fresh or frozen samples that have been stabilized with EDTA or citrate. Erythrocytes are removed in an initial lysis step. After an second lysis step, genomic DNA is bound to a Mini Filter DNA, which can be discarded. RNA is selectively bound to a Mini Filter RNA, washed with two different buffers and eluted. Including lysis, isolated RNA is available in approximately 45 min. The isolated RNA is suitable for a wide range of different molecular biology applications, including RT-PCR.

The GenUP Blood RNA Kit is designed for the use with blood. For other starting material, such as cell-free body fluids (including cerebrospinal fluid, serum, plasma or urine), tissue, stool samples, buffy coat, cultured or isolated cells, swabs, dried blood spots, viruses, fungi, bacteria or parasites, please refer to the GenUP Total RNA Kit (BR07009), GenUP Virus RNA Kit (BR07010) or GenUP Virus DNA/RNA Kit (BR07011).

DESCRIPTION

biotechrabbit[™] GenUP[™] Plant RNA Kit has been developed for quick and easy purification of total RNA from plant materials. After initial homogenization and lysis, genomic DNA is bound to a Mini Filter DNA, which can be discarded. RNA is selectively bound to a Mini Filter RNA, washed with two different buffers and eluted. The purified RNA is ready for use in any demanding molecular biology application, including RT-PCR.

Two lysis buffers, Buffer LYSIS LR and Buffer LYSIS LT, are provided to maximize yield. Most plant material can be processed with Buffer LYSIS LR. In the cases that yield using Buffer LYSIS LR is low, use Buffer LYSIS LT.

GenUP™ BS Virus RNA Kit

Cat. no.	Size	Package content
BR0702303	250 preps	Solutions, Mini Filters
		and Vials
		-
Starting material Swab samples		
Extraction time	Approximately 25 min	
Typical yield	Yield is dependent on sample quality	

FEATURES

- Fast and simple procedure
- High yields of pure RNA
- No DNase treatment, no toxic β-mercaptoethanol

and viral load

APPLICATIONS

Virus RNA isolation from swabs

DESCRIPTION

biotechrabbit[™] GenUP[™] BS Virus RNA Kit has been specially developed for quick and easy isolation of viral RNA from swabs, including nasopharyngeal and oropharyngeal swabs. The unique binding membrane of our high-capacity Mini Filters guaranties high yields. A high concentration of purified RNA can be achieved with flexible elution volumes. The kit includes carrier RNA.

After few initial steps, the viral RNA is bound to a Mini Filter, washed, and then eluted in a separate tube. The purified RNA is ready to be used in all demanding molecular biology applications, including qRT-PCR.

GenUP[™] Virus DNA/RNA Kit

Cat. no.	Size	Package content	
BR0701101	10 preps	Solutions, Mini Filters	
BR0701102	50 preps	and Vials	
BR0701103	250 preps		
		<u> </u>	
Starting material Eukaryotic cells (up to 5 × 10 ⁶),			
	serum, plasma, cell-free body fluids,		
	cell culture supernatants (150 µl), tissue		
	samples, biopsies (up to 20 mg), paraf-		
	fin-embedo	fin-embedded tissues, buccal swabs	
Extraction time	Approximately 25 min		
Typical yield	Yield is hig	hly dependent on sample	
	type		

FEATURES

- Universal kit for simultaneous RNA and DNA isolation from different starting materials
- · Fast and simple procedure, sample specific protocols
- High yields of pure RNA and DNA
- No DNase treatment, no toxic β-mercaptoethanol

APPLICATIONS

- Simultaneous viral DNA and RNA isolation from various sources
- Excellent performance for unknown viruses

DESCRIPTION

biotechrabbit[™] GenUP[™] Virus DNA/RNA Kit has been specially developed for quick and easy isolation of viral RNA and DNA. The kit is especially useful when the origin of the virus is unknown. Viral double-stranded DNA and singlestranded RNA are simultaneously isolated from eukaryotic samples, including plasma, serum, and other body fluids as well as cell cultures, tissues, and buccal swabs.

The unique binding membrane of our high-capacity Mini Filters guaranties high yields. A high concentration of purified nucleic acid can be achieved with flexible elution volumes. The kit includes carrier RNA.

After a few initial procedures, the viral nucleic acids are bound to a Mini Filter, washed and then eluted in a separate tube. The purified nucleic acids are ready to be used in all demanding molecular biology applications, including cDNA synthesis, northern blot analysis, qPCR and RT-PCR.



GenUP[™] FFPE Paraffin Removal Solution

Cat. no.	Size	Package content	
BR0701601	10 ml	GenUP FFPE Paraffin	
BR0701602	50 ml	Removal Solution	
Starting material 2–4 fresh slices from a FFPE sample			
Extraction time	Approximately 45–60 min		
Composition	Single solution formulation for fast		
	and easy DNA extraction		
DNA isolation	Compatible with all common genomic		
	DNA purific	DNA purification kits	

FEATURES

- One-step deparaffinization of FFPE slices
- No toxic or flammable chemicals

APPLICATIONS

• Removal of paraffin from formalin-fixed, paraffin-embedded (FFPE) sample slices

DESCRIPTION

biotechrabbit[™] GenUP[™] FFPE Paraffin Removal Solution is an odor-free, biofriendly and innovative solution to deparaffinize FFPE slices within minutes in a single step. This product replaces conventional procedures that involve toxic and/or flammable solvents. After removal of paraffin from FFPE slices with GenUP FFPE Paraffin Removal Solution, continue with the genomic DNA purification kit of your choice.

Electrophoresis

Stable and ready-to-use, biotechrabbit™ DNA and protein ladders provide dyes, allowing the progress of gel electrophoresis to be monitored. The location and intensity of ladder bands in gels can be used to estimate the size and, for DNA, quantity of other bands in the gel. Nuclease-free DNA Loading Dye is also available.

- Time saving ready to use for convenience
- Reliable high purity and stability

DNA Electrophoresis Ladders and Loading Dye

Cat. no.	Size	Package content	
1 kb DNA Ladder with 6X Loading Dye			
BR0800101	100 lanes	 1 kb DNA Ladder RTU 	
		 6X DNA Loading Dye 	
100 bp DNA Ladder with 6X Loading Dye			
BR0800201	100 lanes	 100 bp DNA Ladder RTU 	
		 6X DNA Loading Dye 	
50 bp DNA Ladder with 6X Loading Dye			
BR0800401	100 lanes	 50 bp DNA Ladder RTU 	
		 6X DNA Loading Dye 	
6X DNA Loading Dye			
BR0800301	5 ml	 6X DNA Loading Dye 	

FEATURES

- · Ready to use DNA ladders ideal for DNA sizing and gel quantification
- Pure and stable retain sharp bands after 6 months storage at room temperature
- Supplied with 6X Loading Dye for sample DNA

APPLICATIONS

· DNA sizing and approximate quantification on agarose gels

DESCRIPTION

biotechrabbit™ DNA electrophoresis ladders are mixtures of exceptionally purified DNA fragments created either by PCR or by digesting proprietary plasmids with restriction enzymes. Ladders are ready to use and suitable not only for DNA sizing but also for approximate DNA quantification in gels. For convenience, ladders have increased intensity reference bands and indicated DNA amount in nanograms for every band.

Every ready-to-use ladder is supplied with the nucleasefree Loading Dye Solution, which ensures optimal migration and quantification of your DNA probes. It includes three electrophoresis tracking dyes (xylene cyanol, bromophenol blue and orange G), allowing the process of the DNA through the gel to be visualized.

Ready-to-use ladder: Range | bands: Reference:

1kb DNA Ladder 250-10,000 bp | 13 1,000 and 3,000 bp

Base Pairs

6,000

4,000

- 3,000 - 2,500

2.000

1.500

- 1.000

- 750

- 500

- 250

DNA Mass

(ng/5µl)

92 34

34 20

92

23

30

45





200 and 500 bp

6X DNA Loading Dye Migration of dyes in 1% TAE agarose gel



(ng/5µl)

DNA Mass




Protein Electrophoresis Ladders

Cat. no.	Size	Package content
TriColor Broad Protein Ladder (3.5–245 kDa)		
BR0900101	100 lanes	 TriColor Broad Protein
		Ladder (3.5–245 kDa)
TriColor Protein Ladder (10–180 kDa)		
BR0900201	100 lanes	TriColor Protein Ladder
		(10–180 kDa)

FEATURES

- Ready-to-use prestained protein ladders covering 3.5–245 kDa and 10–180 kDa ranges
- Pure and stable retain sharp bands after 3 months storage at 4 °C
- Protein bands in three colors with easily recognizable colored reference bands

APPLICATIONS

- Approximate protein sizing on SDS-PAGE and western blots
- Monitoring protein gel electrophoresis and western transfers

DESCRIPTION

biotechrabbit[™] TriColor Broad Protein Ladder facilitates approximate molecular-weight estimation of proteins on denaturing polyacrylamide gels, monitoring protein separation during electrophoresis and verification of western transfer efficiency to membranes (polyvinylidene difluoride, nylon or nitrocellulose). The ladder provides three colored reference bands.

TriColor Broad Protein Ladder is ready for immediate use and is notably stable: sharp bands are produced after storage for 2 weeks at room temperature or 3 months at 4 °C, eliminating the need to thaw before loading.

Ready-to-use ladder:

Range | bands: Reference: TriColor Broad Protein Ladder (3.5–245 kDa) 3.5-245 kDa | 13 25 and 75 kDa

TriColor Protein Ladder (10–180 kDa) 10–180 kDa | 10 25 and 75 kDa



Migration patterns in different electrophoresis conditions. The apparent molecular weight of each protein (kDa) was determined by calibrating against unstained proteins under same conditions.

NOTES



Cell-Free Protein Synthesis

Cell-free protein synthesis overcomes the limitations of traditional cell-based methods. Advantages are ease of use, speed, throughput and the flexible character of the system. Expression conditions can be controlled and easily adapted by adding supplements. It is possible to express toxic proteins and to incorporate labeled amino acids into the protein of interest.

Laborious up- and downstream procedures of traditional *in vivo* methods are avoided as this technique does not require gene transfection or cell culture. By using linear PCR based templates to express the gene of interest, time-consuming cloning procedures are completely avoided.

biotechrabbit[™] expression systems are based on extracts from *E. coli*, wheat germ and insect cells. All components are optimized for highest productivity.

- · From gene to protein within one day
- No transfection, no cell culture
- Expression of toxic proteins
- · Easy site-specific labeling of proteins
- Excellent reproducibility
- High-throughput screening approaches
- Dedicated His-tag protein purification kits

Cell-Free Expression Workflow

In vitro transcription and translation



Protein analysis or functional assay

Continuous Exchange Cell-Free (CECF) Technology

Cell-free protein expression in batch format is highly suitable for screening applications because of the fast and easy setup. However, due to inhibition from byproducts and depletion of substrates, yields using this process are limited. Highest yields are achieved with continuous exchange during cell-free expression (CECF), which allows protein expression to continue for 24 hours. CECF relies on diffusion through a semipermeable membrane between the reaction and the feeding chamber. The reaction chamber contains the machinery for protein expression, while the feeding chamber continuously supplies amino acids, NTPs and energy-regenerating compounds. During the reaction, inhibitory byproducts diffuse out off the reaction chamber. In combination with the optimized high-yield (HY) lysates, synthesis rates using CECF are much higher compared to conventional cell-free expression systems.

Batch







Substrates

Semipermeable membrane

RTS™ 100 E. coli HY Kit

Cat. no.	Size	Package content
BR1400106	6 rxns of 50 µl	Reagents for coupled
BR1400101	24 rxns of 50 µl	transcription/translation
BR1400102	96 rxns of 50 µl	reactions

Proteins expressed	Standard soluble proteins
	Membrane proteins
Reaction time / yield	1–4 h / 20 µg
Reaction mode	Batch
Template	RTS Linear Template Kit Plus
	RTS Linear Template Fab Kit
	RTS 100 E. coli LinTempGen Set, His-tag
	RTS pIX3.0 Vector
	RTS pIVEX <i>E. coli</i> His-tag, 2nd Gen.

FEATURES

- · Compatible with linear templates for fast screening without cloning
- · Convenient reaction setup
- Fast expression of proteins in parallel
- Improved high yield biochemistry for up to 20 µg protein per reaction

APPLICATIONS

• Fast parallel microgram-scale protein expression

DESCRIPTION

The biotechrabbitTM RTSTM 100 *E. coli* HY Kit contains a modified high-yield *E. coli* lysate to drive coupled *in vitro* transcription and translation reactions. The enhanced kit can express up to 20 μ g protein (50 μ l reaction) in 4 hours using a batch mode. The RTS 100 *E. coli* HY kit is compatible with both linear templates and plasmids, making it the product of choice for protein-expression screening and optimization.

Each kit contains everything you need to perform proteinexpression in the delivered microplate or, alternatively, in microtubes.

RTS™ 500 ProteoMaster E. coli HY Kit

Standard caluble proteins

Size	Package content
5 rxns of 1 ml	Reagents for coupled transcription/translation reactions
	Size 5 rxns of 1 ml

Froteins expressed	Stanuaru soluble proteins
	Membrane proteins
Reaction time / yield	6-24 h / 6 mg
Reaction mode	CECF
Template	RTS pIX3.0 Vector
	RTS pIVEX E. coli His-tag, 2nd Gen.

FEATURES

Drataina avaragad

- Unique CECF technology to ensure high yields
- Enable expression of toxic proteins that cannot be produced *in vivo*
- Incorporation of labeled methionine into synthesized proteins

APPLICATIONS

- Optimized medium-scale, cell-free expression of up to 6 mg protein
- Protein synthesis for NMR spectroscopy or X-ray crystallography

DESCRIPTION

The biotechrabbit[™] RTS[™] 500 ProteoMaster *E. coli* HY Kit is designed and optimized for the expression of up to 6 mg of protein from plasmid templates in highly efficient coupled transcription/translation reactions. This is achieved by utilizing a biochemically enhanced high-yield (HY) *E. coli* lysate and the RTS 500 ProteoMaster reaction device.

Due to preparative expression yields, the RTS 500 ProteoMaster *E. coli* HY Kit is suited for the production of protein for NMR spectroscopy or X-ray crystallography. The kit design allows to use amino-acid mixtures of your choice. Methionine (included in the kit separately from other amino acids) can be easily substituted by selenomethionine (not included). Other amino acids can be exchanged by preparing tailored amino-acid mixtures using the RTS Amino Acid Sampler.

The RTS 500 ProteoMaster *E. coli* HY Kit is optimized use in a thermomixer. The continuous agitation of the device guarantees optimal support of the CECF process, which is required for efficient protein expression.



RTS™ 9000 E. coli HY Kit

Cat. no.	Size	Package content
BR1400301	1 rxn of 10 ml	Reagents for coupled transcription/translation reactions
Proteins expr	essed Standar	d soluble proteins

	Membrane proteins
Reaction time / yield	6–24 h / 50 mg
Reaction mode	CECF
Template	RTS pIX3.0 Vector
	RTS pIVEX E. coli His-tag, 2nd Gen.

- FEATURESUnique CECF technology to ensure high yields
- Expression of toxic proteins that cannot be produced in vivo
- · Incorporation of labeled methionine into proteins

APPLICATIONS

• Large-scale, cell-free expression of up to 50 mg protein

DESCRIPTION

One reaction of the biotechrabbit[™] RTS[™] 9000 *E. coli* HY Kit yields ten times more total protein (up to 50 mg) than an RTS 500 reaction. The kit provides a reaction volume ten times higher than the RTS 500 *E. coli* HY Kit, while maintaining a similar productivity per milliliter of lysate.

The RTS 9000 Reaction Device is composed of a reaction compartment and a feeding compartment. The 10 ml reaction compartment, where the coupled transcription/ translation reactions take place, is separated from the 100 ml feeding compartment by a semipermeable membrane (cut-off 10 kDa).

RTS™ 100/500 E. coli Disulfide Kits

Cat. no.	Size	Package content
RTS 100 <i>E. coli</i> Disulfide Kit		
BR1400401	24 rxns	Reagents for coupled transcription/
	of 50 µl	translation reactions
RTS 500 E. coli Disulfide Kit		
BR1400501	5 rxns	Reagents for coupled transcription/
	of 1 ml	translation reactions
Proteins expr	ressed	Standard soluble proteins
		Disulfide-bonded proteins
		(e.g., Fab antibodies)
Reaction time / yield		6–24 h / 80 µg (50 µl), 2.5 mg (1 ml)
Reaction mode		CECF
Template		RTS pIX3.0 Vector
		RTS pIVEX <i>E. coli</i> His-tag, 2nd Ger

FEATURES

- Optimized reaction conditions for the expression of functional disulfide-bonded proteins
- No need for refolding aggregated proteins
- CECF technology for yields ables of up to 2.5 mg

APPLICATIONS

- Expression of active proteins containing multiple disulfide bonds
- Expression of Fab antibodies for rapid functional testing

DESCRIPTION

The biotechrabbit[™] RTS[™] 100/500 *E. coli* Disulfide Kits provide an optimized combination of lysate, reagents, chaperones, and buffer systems to produce maximum yields of active proteins containing multiple disulfide-bonds in a CECF system.

The kit employs several innovative technologies:

- Optimized high-yield (HY) E. coli lysate
- Redox buffer to maintain the system under oxidizing conditions, which enhances disulfide-bond formation
- Disulfide isomerase for rearrangement of disulfide bonds, which enhances the yields of correctly folded proteins

The CECF technology prolongs the reaction time up to 24 hours and, therefore, enhances protein yield. Transcription, translation and disulfide-bond formation take place in the reaction compartment of the device. Substrates and energy components needed for a sustained reaction are continuously supplied via a semipermeable membrane. At the same time, potentially inhibitory interactions with byproducts are diluted by diffusion through the same membrane into the 1ml feeding chamber. Depending on the particular protein, expression of up to several milligrams of protein per reaction can be obtained (e.g., > 2 mg/ml of active urokinase).

RTS[™] 100 Wheat Germ Kit

Cat. no.	Size		Package content
BR1402501	24 rxns o	of	Reagents for coupled
	50 µl		transcription/translation
			reactions
Proteins expre	essed	Stand Com	dard soluble proteins blex and large proteins
Reaction time Reaction mod Template	/ yield e	Mem 3h/5 Batch BTS	brane proteins 5 µg 1 Wheat Germ LinTempGen Sei
		RTS	olVEX Wheat Germ His ₆ -tag

FEATURES

- Efficient expression of up to 5 µg protein per reaction
- Compatible with non-optimized, pro- and eukaryotic T7 promoter-based vectors as well as linear templates
- Excellent homogeneity of large proteins
- No endotoxins, allows direct use in cell-based assays
- Upscaling using high yield CECF wheat germ kits

APPLICATIONS

- Fast parallel synthesis for screening purposes
- Synthesis of mammalian or other proteins for functional testing without the disturbing endogenous protein background that is typical for *in vivo* expression
- Synthesis of proteins dependent on eukaryotic chaperones

DESCRIPTION

The biotechrabbit[™] RTS[™] 100 Wheat Germ Kit is a batch screening system allowing rapid protein expression within 2.5 hours. The kit allows use of a broad range of non-optimized T7 promoter-based expression vectors from prokaryotic and eukaryotic origin in addition to established cell-free wheat germ vectors or linear templates. Using this approach, multiple constructs — including vectors that have not been optimized for wheat germ — can be screened rapidly for expression and functionality in parallel. The procedure saves precious time otherwise required for template generation. The new batch system yields up to 100 µg protein per ml reaction, which is sufficient for many downstream applications.

RTS[™] 100/500 Wheat Germ CECF Kits

Cat. no.	Size	Package content
RTS 100 Wh	eat Germ	CECF Kit
BR1401001	24 rxns	Reagents for coupled transcription/
	of 50 µl	translation reactions
RTS 500 Wheat Germ CECF Kit		
BR1401101	5 rxns	Reagents for coupled transcription/
	of 1 ml	translation reactions
Proteins exp	ressed	Standard soluble proteins
		Complex and large proteins
		Membrane proteins
Reaction time / yield		6–24 h / 50 µg (50 µl), 1 mg (1 ml)
Reaction mode		CECF
Template		RTS Wheat Germ LinTempGen Se
		RTS pIVEX Wheat Germ His6-tag

FEATURES

- Successful expression of eukaryotic proteins, generally without sequence optimization
- Synthesis of up to 1 mg functional eukaryotic protein
- Maximized production of active heat-sensitive proteins, due to an optimum reaction temperature of 24 °C

APPLICATIONS

- Up to milligram-scale protein expression of functional eukaryotic proteins
- Expression of multiple proteins in parallel

DESCRIPTION

The biotechrabbit[™] RTS[™] 100/500 Wheat Germ CECF Kits enable high success rates for expression and solubility, particularly for eukaryotic target proteins. The use of PCRgenerated linear templates greatly improves convenience. The kits allow synthesis of proteins in a wide molecular weight range (up to 220 kDa, as tested), as well as synthesis of truncated protein variants (e.g., for epitope or functional domain mapping). While the RTS 100 Wheat Germ CECF Kit is designed for high-throughput protein expression, the RTS 500 version ensures higher yields.

The system can be used with linear templates as well as mRNA, vectors and plasmid DNA. Depending on the protein, expression yields of more than 1 mg protein per reaction can be obtained within 24 hours in a 1 ml reaction.



RTS[™] 100 Insect Membrane Kit

Cat no	Size	Package content
BR1401501	5 ryps of 50 ul	Reagents for linked
DR1401501		
BR1401502	20 rxns of 50 µl	transcription/translation
		reactions
Proteins expres	ssed S	andard soluble proteins
	С	omplex and large proteins
	G	lycosylated proteins
	Ν	embrane proteins
Reaction time /	yield 4	h/2µg
Reaction mode	B	atch
Template	R	TS Linear Template Kit Plus
	R	TS Linear Template Fab Kit
	R	TS pIX4.0 Insect Vector
	R	TS pIX3.0 Vector
	R	TS pIVEX Wheat Germ Hise
	R	TS DIVEX E coli His-tag 2nd
		TS 100 F calil in Tomp Can
	K	15 IUU E. COII LINTempGenS

APPLICATIONS

- Expression of membrane proteins without the need for additives
- Synthesis of post-translationally modified eukaryotic proteins

RTS[™] Amino Acid Sampler

Cat. no.	Size	Package content
BR1401801	5 rxns of 1 ml	20 L-amino acids (1 vial each)
		5 × DTT (1,4-Dithiothreitol)

FEATURES

• Set of 20 individual amino acids, allowing amino-acid specific protein labeling for structure determination

APPLICATIONS

- For use with RTS 100, RTS 500 and RTS 9000 E. coli kits
- · Ideal for labeling of proteins for NMR and X-ray analyses

FEATURES

- Robust expression of a wide range of proteins (clotting factors, protein kinases, transcription factors)
- Yields of up to 50 µg/ml protein, sufficient for functional studies or screening procedures
- Compatibility with linear templates and plasmids
- Expression screening in small volumes (> 5 µl)

DESCRIPTION

The biotechrabbit[™] RTS[™] 100 Insect Membrane Kit was the first commercially available *in vitro* protein synthesis kit based on cell lysates derived exclusively from insect cell cultures. These homogeneous lysates from a Spodoptera frugiperda cell line enable efficient expression of a broad range of eukaryotic proteins with posttranslational modifications. Now, with an optimized lysate production procedure, this kit is even more efficient.

In addition, the lysates contain functional organelle membrane fractions, whose activity is required for posttranslational modification of eukaryotic proteins, including membrane proteins.

RTS 100 Insect Membrane Kit can be used with linear templates as well as expression vectors.

DESCRIPTION

Using RTS[™] 500 ProteoMaster or RTS 9000 *E. coli* HY Kits, protein yields are sufficient in many cases for subsequent use in NMR spectroscopy. The RTS Amino Acid Sampler provides ready-to-use stock solutions of amino acids which can be labeled and used to replace the amino acids provided in the RTS *E. coli* HY Kits.

RTS[™] DnaK Supplement

Cat. no.	Size	Package content
BR1401601	5 rxns of 1 ml	DnaK Mix (DnaK Supplement)
		Energy Mix (DnaK Supplement)

FEATURES

- *E. coli* chaperones to enhance correct folding and increase solubility, even of large proteins
- Compatible with RTS 100/500/9000 E. coli HY Kits

APPLICATIONS

- · Enhance the expression of soluble proteins
- · Improved synthesis of functionally activity proteins

DESCRIPTION

The Rapid Translation System (RTS) is an "open" cell-free protein expression system, in which reaction conditions can easily be adapted in a protein-specific manner by adding co-factors, ligands, detergents, or chaperones to the reaction mixture.

The biotechrabbit[™] RTS[™] DnaK Supplement includes DnaK (heat shock protein 70, hsp 70), one of the key chaperone systems in *E. coli*, in combination with DnaJ (hsp 40) and GrpE. The RTS DnaK Supplement can be used to increase the amount of correctly folded and/or soluble proteins expressed in the RTS.

DnaK can be used with proteins of any size and serves as an aggregation inhibitor that binds unfolded protein or transfers the protein to other chaperones, such as the GroE or the Hsp90 systems. The RTS DnaK Supplement can be used with the RTS *E. coli* HY platform.

RTS™ GroE Supplement

Cat. no.	Size	Package content
BR1401701	5 rxns of 1 ml	GroEL and GroES
		(GroE Supplement)

FEATURES

- E. coli chaperone mix to enhance correct folding and increase solubility of expressed proteins
- Compatible with RTS 100/500/9000 E. coli HY Kits

APPLICATIONS

Improved synthesis of functionally active proteins

DESCRIPTION

The Rapid Translation System (RTS) is an "open" cell-free protein expression system, in which reaction conditions can easily be adapted in a protein-specific manner by adding co-factors, ligands, detergents, or chaperones to the reaction mixture.

The biotechrabbitTM RTSTM GroE Supplement provides the GroEL and GroES (together referred to as GroE), which are cytoplasmic chaperones of *E. coli* reported to be involve in the de novo folding of 10–30 % of all cytoplasmic proteins. It is not possible to predict that a particular protein is a substrate for GroE, GroEL or GroES.

However, the presence of GroE during cellular or cell-free protein synthesis has been reported to lead to higher amounts of correctly folded and active product in many cases. The RTS GroE Supplement can be used with the RTS *E. coli* HY platform.



Linear Templates and Vectors for Cell-Free Protein Synthesis

biotechrabbit[™] cell-free expression kits can be used with linear templates and expression vectors to produce the proteins of interest. The ability to use linear templates is one of the major advantages of cell-free compared to cell-based expression systems. Linear expression templates can be rapidly generated by PCR, allowing multiple constructs to be screened quickly and easily.

Expression vectors enable protein production to be upscaled. Either the genes of interest are directly inserted into an expression vector or linear templates are generated first by PCR and subsequently cloned. Protein yields achieved are sufficient for extended functional and structural analysis, or production of labeled proteins for NMR spectroscopic or X-ray crystallographic studies.

- Fast PCR-based generation of linear templates
- Rapid expression screening and optimization
- Optimized expression vectors for scaling
 production up to 6 mg per ml
- Flexible positioning of purification tags

No.	Product	Protein tag	Expression kits (see page 61)
	Linear template Kits		
1	RTS Linear Template Kit Plus	His-tag, Strep-tag, Signal peptide	RTS 100 <i>E. coli</i> HY Kit RTS 100 Insect Membrane Kit
2	RTS Wheat Germ Lin TempGen Set, His_6 -tag	His-tag	RTS 100/500 Wheat Germ Kits
3	RTS Linear Template Fab Kit	His-tag, Signal peptide	RTS 100 <i>E. coli</i> HY Kit RTS 100 Insect Membrane Kit
4	RTS 100 <i>E. coli</i> LinTempGen Set, His-tag	His-tag	RTS 100 <i>E. coli</i> HY Kit RTS 100 Insect Membrane Kit
	T7 vectors		
5	RTS pIX4.0 Insect Vector	Tag from cloned gene	RTS 100 Insect Membrane Kit
6	RTS plX3.0 Vector	Tag from linear template	RTS 100/500/9000 <i>E. coli</i> HY Kits RTS 100/500 <i>E. coli</i> Disulfide Kits RTS 100 Insect Membrane Kit
7	RTS pIVEX Wheat Germ ${\rm His}_6{\rm -tag}$ Vector Set	His-tag	RTS 100/500 Wheat Germ Kits
8	RTS pIVEX <i>E. coli</i> His-tag, 2nd Gen. Vector Set	His-tag	RTS 100/500/9000 <i>E. coli</i> HY Kits RTS 100/500 <i>E. coli</i> Disulfide Kits RTS 100 Insect Membrane Kit

RTS[™] Linear Template Kit Plus

Cat. no.	Size	Package content
BR1402401	20 PCRs of 25 µl	PCR reagents for linear
		template generation

Protein tag Expression kits His-tag, Strep-tag, Signal peptide RTS 100 *E. coli* HY Kit RTS 100 Insect Membrane Kit

FEATURES

- Efficient in vitro expression (E. coli and insect kits)
- Primers for affinity tags (His₆-tag /Strep-tag II)
- Signal peptide sequence primers for expression translocation into microsomes (insect systems)

APPLICATIONS

- Fast generation of multiple template constructs
- Accelerated screening of protein expression templates
 and conditions

DESCRIPTION

The RTSTM Linear Template Kit Plus accelerates the generation of multiple templates for parallel protein expression. Constructs are suitable for expression in *E. coli* and insect cell-free lysates. Due to compatible restriction sites in the adapter primers, products can be easily cloned into pIX3.0 vectors for large-scale *in vitro* expression.

The kit contains primers providing regulatory elements required for optimal transcription and translation in cell-free expression systems. Specially designed 5' untranslated regions reduce the formation of secondary structure in the translation initiation region, one of the most common causes of low expression rates. A His-tag or Strep-tag II can be added to either terminus, greatly simplifying protein purification and detection.

A signal sequence enables expression of functional proteins containing disulfide-bonds or membrane proteins in the RTS 100 Insect Membrane kit. A signal peptide can be introduced using the N-terminal primer, which contains a melittin sequence. This enables protein translocation into microsomal vesicles derived from insect cell endoplasmic reticulum.[™]

RTS[™] Wheat Germ LinTempGen Set, His₆-tag

BR1401201 96 PCRs of 50 µl PCR reagents fo	or linear
template generation	ition

His-tag RTS 100 Wheat Germ Kit RTS 100 Wheat Germ CECF Kit RTS 500 Wheat Germ CECF Kit

FEATURES

Protein tag

Expression kits

- · Optimized spacing of eukaryotic regulatory elements
- Time saving, no laborious cloning
- Flexible positioning of His6-tag at the C- or N-terminus
- · Primary PCR product can be cloned into pIVEX vectors

APPLICATIONS

- Rapid generation of linear expression constructs for screening applications
- Linear template generation for optimized His₆-tagged protein expression

DESCRIPTION

The RTS[™] Wheat Germ LinTempGen Set, His6-tag can be used to rapidly generate linear expression constructs that are ready for use with RTS Wheat Germ kits. PCR generated linear templates allow fast protein expression from wild type or modified DNA sequences. Use PCRbased template generation to:

- · Add epitope-tag sequences for antibody detection
- Introduce mutations
- Change codon usage
- Generate truncated proteins

Using linear templates for protein expression avoids the time-consuming step of cloning into expression plasmids and can be used for initial, rapid screening for expression. Gene-specific primers can be used for RTS Wheat Germ and *E. coli* LinTempGen sets enabling easy transfer of products from any expression system to both the RTS Wheat Germ and *E. coli* platforms. For expression scale-up, the linear templates can be cloned into the RTS plVEX Wheat Germ vector and used with RTS 100/500 Wheat Germ CECF Kits.



RTS[™] Linear Template Fab Kit

Cat. no.	Size	Package content
BR1402201	96 PCRs of 25 µl	PCR reagents for linear
		template generation

Protein tag Expression kits His-tag, Signal-peptide RTS 100 *E. coli* HY Kit RTS 100 Insect Membrane Kit

FEATURES

- Introduction of an affinity tag for protein purification
- Signal peptide sequence for translocation into microsomes
- Introduction of a signal peptide sequence for translocation into microsomes
- · Compatibility with E. coli HY and Insect kits

APPLICATIONS

- PCR-based inear templates for protein expression in *E. coli* and insect kits
- Fast template generation for expression of Fab and scFv antibody fragments

DESCRIPTION

The RTS[™] Linear Template Fab Kit is designed for rapid PCR-based generation of templates directly used in the *E. coli* Disulfide and Insect kits allowing functional screening of antibody fragments and disulfide proteins and to facilitate the assignment of biological functions to genes. The templates can be generated with optional affinity tag and/ or signal peptide sequences for protein purification and translocation into microsomes.

The RTS Linear Template Fab Kit generates templates with elements to drive efficient coupled *in vitro* transcription and translation reactions for the synthesis of proteins with intraand intermolecular disulfide bonds. The kit is the product of choice for functional screening of disulfide proteins and optimization of constructs. One kit contains sufficient volumes of primer solutions to perform 96 standard protein expression reactions with reaction volumes of 25 μ l.

RTS™ 100 E. coli LinTempGen Set, His-tag

Cat. no.	Size	Package content
BR1400801	96 PCRs of 50 µl	PCR reagents for linear
		template generation
Protein tag	His-ta	a

Expression kits

His-tag RTS 100 *E. coli* HY Kit RTS 100 Insect Membrane Kit

FEATURES

- · Contains all regulatory elements in optimal spacing
- Option for templates with N- or C-terminal His₆-tag
- · Primary PCR product can be cloned into pIVEX vectors
- For use with RTS 100 *E. coli* HY and insect kits

APPLICATIONS

- Fast gene template generation without cloning
- Generate linear templates for rapid screening and optimization of His₆-tagged protein expression

DESCRIPTION

The RTS[™] 100 *E. coli* LinTempGen Set is designed to rapidly generate linear expression constructs that are ready for use with RTS 100 *E. coli* HY and insect kits. PCR generated linear templates allow fast protein expression from wild type or modified DNA sequences. Use PCR-based template generation to:

- · Add epitope-tag sequences for antibody detection
- Introduce mutations
- Change codon usage
- Generate truncated proteins

Using linear DNA templates for protein expression avoids the time-consuming step of cloning into expression plasmids and can be used for initial, rapid screening for expression. Gene-specific primers can be used for RTS Wheat Germ and *E. coli* LinTempGen sets enabling easy transfer of products from any expression system to both the RTS Wheat Germ and *E. coli* platforms. For expression scale-up, the linear templates can be cloned into RTS pIVEX *E. coli* vectors and used with RTS 100/500/9000 *E. coli* CECF kits.

RTS[™] pIX4.0 Insect Vector

Cat. no.	Size	Package content
BR1400901	1 vector	pIX4.0 Insect Vector

Protein tag Expression kits Tag from cloned gene RTS 100 Insect Membrane Kit

FEATURES

- Regulatory elements for efficient transcription
 and translation
- Multiple cloning site and translational stop codons in all reading frames
- Optimized untranslated regions stabilizing and protecting RNA from degradation by exonucleases
- RNA of defined length via optional plasmid linearization

APPLICATIONS

 Cloning genes for efficient synthesis of proteins using the RTS 100 Insect kits

DESCRIPTION

The RTS[™] plX4.0 Insect Vector is designed for cloning genes into its multiple cloning site. The vector contains efficient untranslated regions specific for the insect expression system, a T7 transcription promoter and additional elements allowing high-yield expression using the RTS 100 Insect Membrane Kit.

The RTS pIX4.0 Insect Vector offers efficient initiation of translation through a T7 promoter element combined with an optimized 3' untranslated region, a T7 terminator, and an optimally positioned linearization site. The expression template is added to the *in vitro* transcription or coupled transcription/translation reactions.

RTS[™] pIX3.0 Vector

Cat. no.	Size	Package content
BR1402701	1 vector	pIX3.0 Vector
Protein tag		Tag from linear template
Expression kit	S	RTS 100 <i>E. coli</i> HY Kit
		RTS 500 ProteoMaster E. coli HY Ki
		RTS 9000 <i>E. coli</i> HY Kit
		RTS 100 <i>E. coli</i> Disulfide Kit
		RTS 500 <i>E. coli</i> Disulfide Kit
		RTS 100 Insect Membrane Kit

FEATURES

- Multiple cloning site compatible with linear templates generated with RTS Linear Template Kit Plus and RTS Linear Template Fab Kit
- High copy origin of replication
- Cell-free protein synthesis scale-up to after expression screening using linear templates
- Free of any potentially interfering regulatory elements ensuring efficient transcription and translation
- · Compatible with RTS E. coli and Insect kits

APPLICATIONS

• Cloning of RTS Linear Templates for expression in RTS *E. coli* and Insect kits

DESCRIPTION

The cloning vector RTSTM pIX3.0 is designed for cloning linear templates generated with RTS Linear Template Plus and Fab kits to subsequently up-scale protein synthesis in *E. coli* batch systems to the CECF format or to use with RTS insect kits. The vector contains a multiple cloning site compatible with that of the linear templates.

The RTS pIX3.0 Vector enables high-level protein expression using the RTS 100/500/9000 *E. coli* HY kits, *E. coli* Disulfide and RTS Insect kits. The kit exhibits the same performance as the RTS pIVEX His-tag, 2nd Gen. Vector Set, which is used for classic cloning.



RTS[™] pIVEX Wheat Germ His₆-tag Vector Set

Cat. no.	Size	Package content
BR1401301	2 vector	pIVEX1.3 WG (His ₆ -tag Vector Set)
		pIVEX1.4 WG (His ₆ -tag Vector Set)
Protein tag Expression ki	its	His-tag RTS 100 Wheat Germ Kit RTS 100 Wheat Germ CECF Kit RTS 500 Wheat Germ CECF Kit

FEATURES

- Stable circular templates ensure high yields in preparative-scale expression reactions
- For construction of C- and N-terminal His₆-tag proteins

APPLICATIONS

• Expression vector for preparative-scale expression in eukaryotic wheat germ lysates

DESCRIPTION

The RTS[™] pIVEX Wheat Germ vectors are designed for high yield cell-free expression of His6-tagged eukaryotic proteins using the RTS Wheat Germ kits.

The two vectors contain all regulatory elements necessary for *in vitro* expression based on a combination of T7 RNA polymerase and wheat germ lysates. They enable the attachment of a His₆-tag in different orientations, at the C- and N-terminus of the protein. Cloning into RTS pIVEX Wheat Germ His₆-tag vectors allows optimal protein expression in RTS Wheat Germ CECF kits to produce more than 1 mg protein per 1 ml synthesis volume in a 24-hour reaction.

RTS™ pIVEX E. coli His-tag, 2nd Gen. Vector Set

Cat. no.	Size	Package content
BR1400701	2 vector	pIVEX2.3d Vector DNA
		2nd Gen. Vector Set
		pIVEX2.4d Vector DNA
		2nd Gen. Vector Set

Protein tag Expression kits His-tag RTS 100 *E. coli* HY Kit RTS 500 ProteoMaster *E. coli* HY RTS 9000 *E. coli* HY Kit RTS 100 *E. coli* Disulfide Kit RTS 500 *E. coli* Disulfide Kit RTS 100 Insect Membrane Kit

FEATURES

- Stable circular templates to ensure high yields in preparative-scale expressions
- For construction of C- and N-terminal His₆-tagged proteins
- Optimal spacing of all regulatory elements
- For use with RTS *E. coli* and Insect kits

APPLICATIONS

Preparative-scale expression of His6-tagged proteins

DESCRIPTION

The RTS[™] pIVEX *E. coli* vectors are designed for high-yield cell-free expression of His₆-tagged proteins using a broad range of RTS expression kits.

The two vectors contain all regulatory elements necessary for *in vitro* expression based on a combination of T7 RNA polymerase and RTS expression lysates. They enable the attachment of a His₆-tag in different orientations at the C- and N-terminus of the protein. Cloning into RTS pIVEX His₆-tag vectors allows optimal protein expression with high yields.

High Affinity His-tag Protein Binding

Modern proteomics relies on reliable and sensitive detection and labeling of tagged proteins, such as those bearing the 6xHis-tag. However, the low affinity of currently available chelators used for protein binding limits the sensitivity of many applications for selective, site-specific conjugation of proteins. To address this issue, biotechrabbit offers a superior His-tag protein ligand for the reversible binding of proteins and cell surfaces: Tris-NTA.

With a binding affinity that is approximately four orders of magnitude higher than monovalent metal ion chelators (such as nitrilotriacetic acid and iminodiacetic acid), Tris-NTA increases the sensitivity in western hybridization, biosensor analysis, functional protein analysis, protein purification and many other applications. 10,000 times higher binding affinity than

conventional chelators

- Reversible binding of biomolecules
- Great flexibility in functional assays



Lata et al., J. Am. Chem. Soc., 2005, 127, 10205–10215

Tris-NTA Amine / Tris-NTA Biotin

Cat. no.	Size	Package content
Tris-NTA Am	ine	
BR1001101	100 µg	Tris-NTA Amine, 1 mg/ml
BR1001102	1 mg	
Tris-NTA Biotin		
BR1001201	100 µg	Tris-NTA Biotin, 1 mg/ml
BR1001202	1 mg	

FEATURES

- A complex of three Ni-NTA groups ensures high-affinity binding of His-tags
- Binding affinity is approximately four orders of
 magnitude higher than monovalent metal ion chelators
- · Protein binding is stoichiometric

APPLICATIONS

- Reversible labeling of proteins or cell surfaces
- Detection and analysis of target molecules
- · Immobilization of proteins, lipids and cells on surfaces
- · Purification and sample preparation of proteins
- Coupling with microscopic or spectroscopic probes



DESCRIPTION

His-tags are one of the most commonly used tags for protein expression analysis. Conventional metal ion chelators, such as nitrilotriacetic acid (NTA) and iminodiacetic acid (IDA), bind His-tags with low affinities in the range of 10 μ M. The biotechrabbit Tris-NTA complexes three NTA groups that together bind a 6xHis-tag with an affinity that is four orders of magnitude higher (1 nM) than is possible with conventional chelators (10 μ M). The binding of His-tags is stoichiometric and single-molecule detection is possible. Binding is reversible: bound His-tags can be released with imidazole or ethylenediamine-tetraacetic acid (EDTA).

biotechrabbit[™] Tris-NTA is available with a free amine group or conjugated to biotin. It can be used in a large range of applications, including protein detection and labeling, coupling proteins, lipids or cells to surfaces, protein purification and reversible protein modification.





NULS

NOTES



Legal Disclaimer and product use limitation

biotechrabbit GmbH makes every effort to include accurate and up-to-date information within this publication; however it is possible that omissions or errors might have occurred. Changes of publication can be made at any time without notice. All mentioned trademarks are protected by law. The purchase of biotechrabbit product does not include a license to perform any patented applications; therefore it is the sole responsibility of users to determine whether they may be required to engage a license agreement depending upon the particular application in which the products are used. biotechrabbit products are developed, manufactured, and sold for *in vitro* use and laboratory use only. They are not suitable for administration to humans or animals. The products are to be used only by trained laboratory personnel who are familiar with potential hazards and in laboratories equipped to perform life science research. biotechrabbit products may contain chemicals which are harmful if misused. Due care should be taken with all products to avoid direct contact with components that are potentially harmful.

Tris-NTA Amine / Biotin

The use of this product for research purposes is covered by a license from QIAGEN Group. No rights to use this product to perform or offer diagnostic, companion diagnostic, commercial testing or other commercial services for money or money's worth are granted by the supply of this product expressly, by implication or by estoppel. Should you wish to use this product for any other purpose not covered by this license, please contact QIAGEN Corporate Business Development at bd@qiagen.com.

Trademarks

AllScript[™], AmPelify[™], AnyScript[™], ApStarTaq[™], biotechrabbit[™], CAPITAL[™], DirectUP[™], ExactiFi[™], FastUP[™], GeneUP[™], RevertUP[™], RTS[™], StranDisplace[™], UniFi[™], UPstart[™], UPzol[™], YourTaq[™] (biotechrabbit GmbH); Agilent[®], Bioanalyzer[®] (Agilent Technologies, Inc.); ROX[™] (Applera Corporation or its subsidiaries); Beacon[®], StepOne[™], Viia[®] (Life Technologies); Bio-Rad[®], CFX384[™], CFX96[™], Chromo4[™], iCycler[®], iq[™], MiniOpticon[™], MyiQ[™], Opticon[™] (Bio-Rad Laboratories, Inc.); Cepheid[®], Smart-Cycler[®] (Cepheid Corp.); Coomassie[®] (Imperial Chemical Industries); Eco[™], Illumina[®] (Illumina Inc.); Eppendorf[®], Mastercycler[®], REALPLEX[®] (Eppendorf AG); LightCycler[®], TaqMan[®] (Roche Group); Mx3000P[®], Mx3005P[®], Mx4000[®], Stratagene[®] (Stratagene); QIAGEN[®], Rotor-Gene[®] (QIAGEN GmbH); Quantica[®] (Jonathan Redfern); Scorpion[®] (DxS Limited); Takara[®], Cycler Dice[™] (Takara Shuzo Co. Ltd.); Techne[®] (Techne Corporation).

www.biotechrabbit.com © October 2024 biotechrabbit GmbH, all rights reserved. PTM-1013-001 valid from 01.10.2024



biotechrabbit GmbH Volmerstr. 9 12489 Berlin, Germany phone: +49 30 555 78 21 0 info@biotechrabbit.com Managing Director: Dr. Bernd Haase Commercial Register: Charlottenburg HRB 232169 B VAT No/USt-Id: DE 279969868

Bank account Deutsche Bank Berlin BIC/SWIFT: DEUTDEDBBER IBAN: DE22100700240062889102